Operating Instructions

F900

Fluorescence Spectrometer Software

Issue 1, March 2004



Edinburgh Instruments Ltd., 2 Bain Square, Kirkton Campus, Livingston, EH54 7DQ. UK T: +44 (0)1506 425300 - F: +44 (0)1506 425320 - E: eai.service@edinst.com - W: www.edinst.com

© Edinburgh Instruments Limited, 2004

All rights reserved. No part of this document may be reproduced or transmitted in any form or by any means without prior written permission of the copyright holder.

The product described in this manual is subject to continuous development and, while every effort has been taken to make sure that the information given is correct, Edinburgh Instruments Limited cannot accept any liability for errors and omissions or their consequences

Table of Contents

1	Introduction					
	1.1 Steady State Fluorimeter (FS920)					
	1.2	Fluorescence Lifetime Spectrometer (FL920)	7			
	1.3	Additional Computer Controlled Options	8			
	1.4	F900 Software Overview	9			
2	Soft	ware Installation	10			
	2.1	Installation of System Disk	10			
	2.2	Installation of Auxiliary Disk	10			
3	Hard	Iware Configuration	11			
	3.1	CD Controller	12			
	3.2	Detectors	12			
	3.3	Lamps	14			
	3.4	Lifetime Cards	16			
	3.5	Miscellaneous	17			
		3.5.1 Multiple Sample Holder	18			
		3.5.2 Oscilloscope	19			
		3.5.3 Cryostat	20			
	36	3.5.4 IIIS Monochromator Tab	20			
	5.0	3.6.1 Stepper Details	23			
		3.6.2 Gratings	24			
		3.6.3 Wavelength Increments	25			
		3.6.4 Ports	25			
	37	3.6.5 Controlled Sills Polarisers	20 27			
4	5.7		21			
4	Son		29			
	4.1	Main Menu and Shortcut Tooldar Status Bar	29			
F	4.2	viring Spectral Data	22			
5	Acq	Anny Spectral Data				
	5.1	Setup Menu: Signal Rate				
			33			
		5.1.2 Using the Signal Rate Screen	33 34			
	5.2	5.1.2 Using the Signal Rate Screen Scan measurements	33 34 34			
	5.2	5.1.2 Using the Signal Rate Screen Scan measurements 5.2.1 Emission Scan	33 34 34 35			
	5.2	 5.1.2 Using the Signal Rate Screen Scan measurements 5.2.1 Emission Scan 5.2.2 Excitation Scan 	33 34 34 35 37			
	5.2	 5.1.2 Using the Signal Rate Screen Scan measurements 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 	33 34 34 35 37 39			
	5.2	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan 	33 34 35 37 39 41 44			
	5.2	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 	33 34 34 35 37 39 41 44 46			
	5.2	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 	33 34 34 35 37 39 41 44 46 46			
	5.2	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 5.3.2 Excitation Correction Scans 	33 34 35 37 39 41 44 46 46 48			
	5.2 5.3 5.4	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 5.3.2 Excitation Correction Scans Map measurements 	33 34 34 35 37 39 41 44 46 46 48 49			
	5.2 5.3 5.4	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 5.3.2 Excitation Correction Scans 5.3.1 Emission Correction Scans 5.3.2 Excitation Correction Scans 5.3.1 Emission Map 5.4.1 Emission Map 	33 34 34 35 37 39 41 44 46 46 48 49 49			
	5.2 5.3 5.4	 5.1.2 Using the Signal Rate Screen Scan measurements 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 5.3.2 Excitation Correction Scans 5.3.2 Excitation Correction Scans 5.4.1 Emission Map 5.4.2 Synchronous Map 5.4.3 Temperature Map 	33 34 35 37 39 41 44 46 46 48 49 49 51 53			
	5.2 5.3 5.4 5.5	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 5.3.2 Excitation Correction Scans Map measurements 5.4.1 Emission Map 5.4.2 Synchronous Map 5.4.3 Temperature Map Anisotropy measurements (performed using optional polarisers) 	33 34 35 37 39 41 44 46 46 48 49 49 51 53 54			
	5.2 5.3 5.4 5.5	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 5.3.2 Excitation Correction Scans 5.3.2 Excitation Correction Scans 5.4.1 Emission Map 5.4.2 Synchronous Map 5.4.3 Temperature Map Anisotropy measurements (performed using optional polarisers) 5.5.1 Emission Anisotropy Scan 	33 34 35 37 39 41 44 46 46 48 49 49 51 53 54 55			
	5.25.35.45.5	 5.1.2 Using the Signal Rate Screen 5.1.2 Using the Signal Rate Screen Scan measurements 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scans 5.3.2 Excitation Correction Scans 5.3.2 Excitation Correction Scans Map measurements 5.4.1 Emission Map 5.4.2 Synchronous Map 5.4.3 Temperature Map Anisotropy measurements (performed using optional polarisers) 5.5.1 Emission Anisotropy Scan 	33 34 35 37 39 41 44 46 46 48 49 49 51 53 54 55 56			
6	5.25.35.45.5Acq	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 5.3.2 Excitation Correction Scans Map measurements 5.4.1 Emission Map 5.4.2 Synchronous Map 5.4.3 Temperature Map Anisotropy measurements (performed using optional polarisers) 5.5.1 Emission Anisotropy Scan 5.5.2 Excitation Anisotropy Scan iring Lifetime Data 	33 34 35 37 39 41 44 46 46 48 49 49 51 53 54 55 56 57			
6	 5.2 5.3 5.4 5.5 Acq 6.1 	 5.1.2 Using the Signal Rate Screen 5.2.1 Emission Scan 5.2.2 Excitation Scan 5.2.3 Synchronous Scan 5.2.4 Kinetic Scan 5.2.5 Cryostat Temperature Scan Correction Scans 5.3.1 Emission Correction Scan 5.3.2 Excitation Correction Scans Map measurements 5.4.1 Emission Map 5.4.2 Synchronous Map 5.4.3 Temperature Map Anisotropy measurements (performed using optional polarisers) 5.5.1 Emission Anisotropy Scan 5.5.2 Excitation Anisotropy Scan 5.5.2 Excitation Anisotropy Scan 	33 34 35 37 39 41 44 46 46 48 49 49 51 53 54 55 56 57 57			

	6.3 Multiple	63
	6.3.1 Sample Set Up Button	64
	6.3.2 Time Set Up Button	65
	6.4 IRES	67
	6.4.2 Time Set Un Button	60 69
	6.5 Anisotropy Lifetime	71
	6.5.1 Sample Set Up Button	73
	6.5.2 Time Set Up Button	74
	6.6 Temperature Lifetime	76
	6.6.1 Sample Set Up Button	77
_	6.6.2 Time Set Up Button	78
7	Viewing Data	81
	7.1 As 2D plot	81
	7.2 As 3D plot	83
	7.3 As Text	86
	7.4 As Contour Plot	86
8	Data Analysis	89
	8.1 Combine (Lifetime and Spectral Data)	89
	8.2 Scale (Lifetime and Spectral Data)	92
	8.3 Normalise (Lifetime and Spectral Data)	92
	8.4 Subtract Baseline (Lifetime and Spectral Data)	94
	8.5 Crop Range (Lifetime and Spectral Data)	94
	8.6 Smooth (Spectral Data)	95
	8.7 Differentiate (Spectral Data)	95
	8.8 Integrate	96
	8.9 Reverse	97
	8.10 Peak Search (Spectral Data)	98
	8.11 Correction (Spectral Data)	99
9	Windows-related Menus	100
	9.1 File Menu	100
	9.2 Edit Menu	104
	9.3 View Menu	104
	9.3.1 Lamp setup	105
	9.3.2 Signal Rate	106
	9.3.3 As 2D Plot, As Contour, As 3D Plot, As Text	106
	9.3.4 Flot options 9.3.5 Zoom In	106
	9.3.6 Cursor	106
	9.3.7 Grid	107
	9.3.8 Peaks	107
	9.3.9 Normalise	107
	9.3.10 Status bai 9.3.11 Tool bar	107
	9.4 Options Menu	107
	9.4.1 Plot defaults	107
	9.4.2 Correction files	107
	9.4.3 Windows options	108
	9.4.4 Hardware configuration	109
	9.4.6 Drivers and cards	110
	9.5 Window Menu	110
	9.6 Help Menu	112
10	Warranty	113
	-	

11 CE Declaration of Conformity

115

1 Introduction

The F900 software is used to control the fluorescent spectrometers manufactured by Edinburgh Instruments Ltd and, in addition, acquire, process and analyse the data produced by them. The instruments are the FS920 Steady State Fluorimeter, the FL920 Fluorescence Lifetime Spectrometer, and the FLS920 Combined Steady State and Lifetime Spectrometer. An overview of steady-state and lifetime options will now be given, together with descriptions of the optional attachments for them.

1.1 Steady State Fluorimeter (FS920)

The FS920 steady state fluorimeter has a modular design. The key components are the xenon lamp, excitation monochromator, sample chamber, emission monochromator, and these individual components are joined together to give a robust opto-mechanical assembly mounted to a platform. Figure 1.1 shows the FS920, comprising xenon lamp (Xe900), excitation and emission monochromators (M300), sample chamber (UC920) with integrated control electronics (CD920/CD930), and detector (S900).

The Xenon lamp produces a broad spectrum of high intensity radiation through the UV-Vis-IR range. The arc from this lamp is imaged onto the entrance slit of the excitation monochromator, which selects a specific excitation wavelength with a specific bandwidth (Figure 1.2).



Figure 1.1: FS920 steady-state fluorimeter configuration.



Figure 1.2: Schematic of the FS920 steady-state spectrometer configuration.

The radiation of a specific bandwidth and spectral resolution is passed from the excitation monochromator through a beam-splitter (where a small percentage is reflected onto a reference detector), an iris (so that the intensity can be controlled) and is then focused into the sample. A portion of any luminescence from the sample is collected by a lens system and focused onto the emission monochromator. The emission monochromator selects a wavelength and spectral bandwidth of this emission and the light is passed into a photomultiplier tube (PMT).

The control electronics are located beneath the sample chamber. Only the xenon lamp power supply and an optional detector cooler power supply are separate from the FS920 spectrometer assembly. This set-up maintains the modularity and flexibility of the opto-mechanical design while minimising the number of external cables. The communication to a separate, dedicated computer is made via an RS232 port. All signal processing electronics are located in the computer. Data acquisition is achieved via a photon counting scalar card (PCS900) plugged into a PCI slot of the computer.

1.2 Fluorescence Lifetime Spectrometer (FL920)

The Fluorescence Lifetime Spectrometer is based on the technique of Time Correlated Single Photon Counting (TCSPC). A typical configuration (Figure 1.3) would include a nanosecond flashlamp (connected to the second port of the excitation monochromator) and the TCSPC data acquisition card, TCC900, which is installed in a PCI slot in the computer. Alternatively, pulsed LED's and pulsed diode lasers can be used as light sources. The light path follows the same route as the FS920 fluorimeter.

Both the FS9290 and the FL920 instruments can be configured as a phosphorescence lifetime spectrometer by using a microsecond lamp, coupled to the first entrance port of the excitation monochromator (also shown in Figure 1.3). The PCS900 computer plug-in card can be set up to operate in multichannel scaling (MCS) mode, a mode acquiring single photons in a time resolved manner.



Figure 1.3: FL920 fluorescence lifetime spectrometer configuration.

1.3 Additional Computer Controlled Options

Along with the standard features of the above instruments, a number of optional components and configurations can be added and controlled by the F900 software.

- Light Sources: Additional light sources can be coupled either directly to the sample chamber or onto the entrance of the excitation monochromator. The F900 software can then control the excitation monochromator, allowing computer control of the excitation source by means of up to two computer controlled swing mirrors. Optional light sources include the nanosecond flashlamp, pulsed LED's or diode lasers, pulsed microsecond flashlamp or high power Q-switched lasers.
- **Detectors:** The monochromators incorporated into the FS920 can be fitted with up to two exit slits. A computer controlled swing mirror enables the user to select the emission port of choice. As such, additional detectors may be fitted to the emission monochromator to increase the detection spectral range. There are currently PMT detectors available that span the wavelength range 185 870 nm. In addition germanium PIN diode detectors can be used with a responsivity from 800 1700 nm or InGaAs PIN photodiodes with responsivity from 900 –2500 nm. Up to 3 diffraction gratings can be mounted on the computer controlled grating turret in the monochromator. The grating is selected to optimise the performance of the monochromator over the wavelength of responsivity of the selected detector. In addition, the PCS900 photon counting card has five signal input channels, each input selected from the F900 software.
- Automated Polarisers: High throughput UV transmitting Glan Thompson polarising prisms, with a 14 mm diameter clear aperture, can be provided to allow anisotropy and other polarisation effects to be studied. The prisms are held in computer controlled mounts allowing user defined or preset polarisation angles.
- **Cryostat Systems:** Oxford Instruments liquid nitrogen or helium cryostats with ITC controllers are used when low temperature measurements are required. The F900 software communicates fully with the cryostat controller allowing temperature profiles and temperature dependant spectral maps to be acquired under computer control.
- **Sample Positioner:** A three position sample holder can be coupled to the sample chamber. This can be controlled by the F900 software and allows different set-ups to be configured for each sample position.

T-Geometry: The system can be configured as a T-Geometry with two emission arms allowing two monochromators and up to four detectors to be permanently fitted to the system.

1.4 F900 Software Overview

The F900 software allows the user to control almost all aspects of the hardware, to set-up measurements, to make measurements, to perform data analysis and process this data. It is a Windows-based operating system and the format of the software is such that functional control can be either from the main menu which can be pulled down to open sub menus or through the shortcut toolbar (Figure 4.1) which opens the sub menus. Each of the highlighted sub menus can open dialog boxes to allow the user to enter data. Sub menus are not active (not highlighted) if the options contained in those sub-menus are not applicable, e.g. the corresponding hardware is not present. Another function of the software is the status bar; this is located at the bottom of the F900 window and displays information about the current setup relevant to the measurement being carried out.

2 Software Installation

The F900 software is the operating software for the FL920, FS920 and the FLS920 spectrometers. The software is compatible with Windows 2000 and it is supplied on two CD-ROMs. The first one contains all the system software and should start automatically when the disk is inserted into the drive. The second CD-ROM contains optional software that has been individually configured for the specific system.

The software has a protection device that needs to be installed in parallel port LPT1 before starting the software.

2.1 Installation of System Disk

To install the F900 software, please follow these instructions:

- 1. Place the F900 installation disk into the CD-ROM drive of the computer. The installation programme should start automatically. If it does not then follow the following procedure:
 - a. On the Windows Taskbar press 'Start' then 'Run'. In this dialog box press 'Browse' and select the CD-ROM drive.
 - b. In the root directory, select the file 'Setup.exe'. Press 'Open' to return to the 'Run' dialog box and then press 'OK'.
- 2. Follow the instructions to install the F900 software package.
- 3. Once the installation has completed, the auxiliary disk for the F900 software can be run.

2.2 Installation of Auxiliary Disk

To install the utilities in the Auxiliary Disk, Please follow the following procedure:

- 1. Place the F900 auxiliary disk into the CD-ROM drive of the computer. The installation programme should start automatically. If it does not then follow the following procedure:
 - a. On the Windows Taskbar press 'Start' then 'Run'. In this dialog box press 'Browse' and select the CD-ROM drive.
 - b. In the root directory, select the file 'Setup.exe'. Press 'Open' to return to the 'Run' dialog box and then press 'OK'.
- 2. In the **Select Features** menu, choose the items to be installed. The options are:
 - a. **Install hardware configuration:** This will load an initial hardware configuration, tailored for the specific system, into Windows registry. Full details of hardware configuration are given in Chapter 3 of this manual.
 - b. **Install correction files:** Checking this option will copy correction files for the instrumentation installed with the system hardware.
 - c. **Install manuals:** Selection of this option will install the manuals relevant to the particular system as portable document format (pdf) files in the EI directory.
 - d. **Install El wallpaper:** Acceptance of this option will install a utility that displays contact information for Edinburgh Instruments.

3 Hardware Configuration

Prior to running a spectrometer with the L900 software, the availability and setting of the spectrometer components have to be made known to the software. This is done by calling the **Hardware Configuration** utility, which loads the spectrometer settings into system Registry for use by the F900 software. This utility can be called from two locations, which are:

- a) The Cd920Test programme that is used in initial configuration of the system and for component testing. This route is only normally used when the system is being initially configured.
- b) The F900 software itself. The hardware configuration utility can be called by selecting Hardware Configuration from the Options menu. This route is normally used when the hardware setup is to be modified for a different type of measurement.

Hardware Device Configuration 🛛 🛛 🗙						
System Type C L Geometry	System Type C L Geometry T Geometry					
Detectors Misc. L	Detectors PC Interface Lifetime Cards Misc. Lamps Monochromators Polarise			me Cards Polarisers		
Excitation Mono Use Grating TM300 I: 200.00 - 900.00 nm					Details	
Emission 1 Mono Use Grating TM300 I: 200.00 - 900.00 nm					Details	
Emission 2 Mono Use Grating TM300 ① 1: 200.00 - 900.00 nm						
Root Directory C:\Program Files\EI\F900						
			OK		Cancel	

Figure 3.1: Opening screen for the Hardware Configuration utility.

When opened, the utility will look similar to the one shown in Figure 3.1. The major components of this screen are the **System Type**, which specifies whether the spectrometer is configured in an **L Geometry** or a **T Geometry**; a set of tabbed items that are groupings for the hardware components (these will be described in more detail below); notification of the root directory for the F900 software; an **OK** button to accept hardware changes that have been made and to write them to system registry; a **Cancel** button to abandon any changes and to exit the utility. The groupings on the tabbed items will now be described in more detail.

3.1 CD Controller



Figure 3.2: Controller tab.

The tab for the CD920/CD930 controller is shown in Figure 3.2. There are two settings in this tab and these are:

Present: Checking this box indicates to the software that the controller is present in the system. This is normally true and it is only for unusual geometries, such as the TL900 lifetime spectrometer, where there is no controller present.

COM Port: This gives the user the choice of RS232 port on the host computer that the controller should be connected . This is normally COM1.

3.2 Detectors

The **Detectors** tab is used to configure the nature of the detectors used to monitor the light output from the Excitation monochromator and to record the light emanating from the emission monochromator(s). An example of this tab is shown in Figure 3.3. It is possible for an emission arm to have more than one detector fitted. This is because the monochromator has two output ports, selectable by a movable mirror.

To inform the software of the presence of a detector, ensure that the appropriate check box has been ticked. For the Reference Detector this is the **Present** check box and for the emission arm detectors these are the **Detector 1** and **Detector 2** check boxes.

Misc.	Lamps	Monochroma	tors	Polarisers				
Detectors	CD Co	ontroller	Li	fetime Cards				
Excitation Refe	Excitation Reference Detector							
Present				Details				
Emission Arm 1								
Detector 1				Details				
Detector 2				Details				
Emission Arm 2								
Detector 1				Details				
Detector 2				Details				

Figure 3.3: Detectors tab.

To enter the details for the Reference Detector press the **Details** button in the **Excitation Reference Detector** group box. This will open a dialog box similar to the one shown in Figure 3.4. The entries that can be made for this box are:

Description: This is a user-chosen name for the reference detector.

Connected to: This is the data acquisition card that is connected to the detector output. The options are the PCS-900 card and the TCC-900 card.

Chan: This is the channel port on the card that is connected to the reference detector.

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

Reference De	etector 🗙
Description:	Ref
Mono Port:	V
Connected to:	PCS-900 💌 Chan: 4 🗼
Default	OK Cancel

Figure 3.4: Example of settings for an excitation reference detector.

To enter the details for an Emission Detector press the **Details** button corresponding to chosen detector's check box. This will open a dialog box similar to the one shown in Figure 3.5. The entries that can be made for this box are:

Description: This is a user-chosen name for the reference detector.

Mono Port: This is the output port of the monochromator to which the detector is attached. The options are Port A and Port B.

Connected to: This is the data acquisition card that is connected to the detector output. The options are the PCS-900 card and the TCC-900 card.

Chan: This is the channel port on the card that is connected to the reference detector.

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

Emission1 De	etector1 🛛 🗙
Description:	R955
Mono Port:	Port B
Connected to:	PCS-900 💌 Chan: 0 🍝
Default	OK Cancel

Figure 3.5: Example of settings for an emission detector.

3.3 Lamps

Detector	s	CD Controller	L L	ifetime Cards
Misc.	Lamps	Monoc	hromators	Polarisers
Nanosecon	d Flashlamp		D	etails
Microsecon	d Flashlamp			1
Present			De	etails
Steady State	e Lamp			1
Present			De	etails

Figure 3.6: Lamps tab.

The Lamps tab is used to configure the light sources that are used to illuminate experimental samples. These lamps are normally fitted to the input ports of the excitation monochromator. An external swing mirror can be

fitted to one of these ports so it is possible to have three alternative light sources fitted to the spectrometer system. If a light source is fitted then the corresponding check box should be ticked.

Nanosecond L	amp 🛛 🔀	
		1
Description:	nF900	
Mono Port:	Port A	
Connected to:	TCC-900 💌 Chan: 0 🍝	
Default	OK Cancel	

Figure 3.7: Example of settings for a nanosecond flashlamp.

To enter the details for a Nanosecond flashlamp press the **Details** button corresponding to the nanosecond flashlamp check box. This will open a dialog box similar to the one shown in Figure 3.7. The entries that can be made for this box are:

Description: This is a user-chosen name for the reference detector.Mono Port: This is the output port of the monochromator to which the detector is attached. The options are Port A and Port B.

Connected to: This is the data acquisition card that is connected to the detector output. The options are the PCS-900 card and the TCC-900 card.

Chan: This is the channel port on the card that is connected to the reference detector.

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

Microsecond Lar	np 🔀
Description	LuE900
Mono Port	Port P
Trigger Delau	
Lamp Latency	
Default	OK Cancel

Figure 3.8: Example of settings for a microsecond flashlamp.

To enter the details for a Microsecond flashlamp press the **Details** button corresponding to the microsecond flashlamp check box. This will open a dialog box similar to the one shown in Figure 3.7. The entries that can be made for this box are:

- **Description**: This is a user-chosen name for the reference detector.
- **Mono Port:** This is the output port of the monochromator to which the detector is attached. The options are Port A and Port B.
- **TriggerDelay:** This is the delay between the trigger pulse to the acquisition card and the one to the flashlamp trigger pulse.
- Lamp Latency: This value is used to delay the time that the acquisition card starts measuring. It compensates for delays between a lamp (or laser) being triggered and its output being generated.

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

Xenon Lamp	×
Description	Xe
Mono Port	Port B
Shutter Device	26
Default	OK Cancel

Figure 3.9: Example of settings for an external lamp.

To enter the details for a steady-state (Xenon) flashlamp press the **Details** button corresponding to the steadystate flashlamp check box. This will open a dialog box similar to the one shown in Figure 3.9. The entries that can be made for this box are:

Description: This is a user-chosen name for the reference detector.

Mono Port: This is the output port of the monochromator to which the detector is attached. The options are Port A and Port B.

Shutter Device: This is the device number used for an external shutter for the steady-state lamp.

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

3.4 Lifetime Cards

Misc.	Lamps	Monochromators		Polarisers
Detectors		CD Controller Li		ifetime Cards
For Lifetime	measuremer	nts use:		
C Norland	I N5000	<u>DLL</u> NONE	<u>Card</u> NONE	
EALTC	● EAI TCC/PCS		DETECTED)

Figure 3.10: Lifetime cards tab.

The **Lifetime Cards** tab displays the availability of the cards that are used for lifetime measurements and if there is more than one, allows the user to choose the card to be used.

3.5 Miscellaneous

Detectors Misc. Lar		CD Controller		Lifetime Cards		
		ps	Monochromate		tors	Polarisers
Multiple Sar Present	nple Hold	der etails etails		Oscillos Pres Oxford 0	cope sent Cryostat sent	Details Details
Temperature	e Sensor					

Figure 3.11: Miscellaneous tab.

The **Miscellaneous** tab is used for settings that do not belong naturally to any of the other tabs. The presence of these attachments is made known to the software by checking the appropriate **Present** box and detailed information for each item can be entered by clicking the corresponding **Details** button.

3.5.1 Multiple Sample Holder

ample Holde	r Configur	ation 🔉
	Description	Default
High Level Info		Low Level Info
Range Low	1	Step Multiplier 9000
Range High	3	Step Zero 0
Initial Position	1	Step Mode Full 💌
Resolution	1	Step Sequence 2 Ph 💌
		Step Freq (ms) 0.99
Default		OK Cancel

Figure 3.12: Example of sample holder settings.

If a multiple sample holder has been installed, then its configuration can be entered on a form similar to the one shown in Figure 3.12. The **High Level Info** grouping refers to the sample holder positions and the **Low Level Grouping** refers to the stepper motor that positions the sample holder. Normally the multiple sample holder has three positions. The low-level settings are not normally changed in the field. Details of the settings are:

- **Description:** This value is entered by the user to put a name to the configuration setup. It is only for identification purposes and has no effect on system operation.
- **Range Low:** This is the lowest available position of the sample holder.

Range High: This is the highest available position of the sample.

- **Initial Position:** This is the position that the sample holder will initially move to when the software is started.
- **Resolution:** This is the increment that is added when the sample holder position is changed. Since the sample holder can only have an integral number of positions this value should always be set to 1.

The **Low Level Info** settings should not be altered from their factory settings, without exlicit instruction from the Edinburgh Instruments service department. The factory settings are:

Step Multiplier: 9000 Step Zero: 0 Step Mode: Full Step Sequence: 2 Ph Step Freq: 0.99

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

3.5.2 Oscilloscope

Oscilloscope Settings	×
Setup	COM Port
Interface: Ethernet	Scope Channel: 1 only
ScopeType: TDS 3012	Data Points: 10000 💌
Default	OK Cancel

Figure 3.13: Example of oscilloscope settings.

If an oscilloscope has been installed with the system, then its configuration can be setup with the form shown in Figure 3.13. The details that can be entered are:

- Interface: This setting is used to select the communications interface between the oscilloscope and the computer. The currently available options are Ethernet and Serial (RS-232). If serial communications have been chosen, the **Setup COM Port** button should be pressed to select the communications values (details below).
- **Scope Channel:** This setting chooses the oscilloscope channel that data is to be acquired from. The options are channel 1 only, channel 2 only, or both channel 1 and channel 2.
- Scope Type: This is the type of oscilloscope that is attached to the system. The currently available options are the Tektronix TDS3012 and Tektronix TDS3052.
- **Data Points**: This is the number of data points that are to be acquired from the oscilloscope. The oscilloscope itself will transfer 10000 data points, but this can be reduced by data compression (averaging of neighbouring data points) to 5000 (1:2), 3333 (1:3), 2500 (1:4), 2000 (1:5), 1000 (1:10), or 500 (1:20). Note that is the time scale on the oscilloscope is less than 200 ns/div then it will always transfer 10000 data points over a time range of 2 μs.

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

If serial communications are to be used, then the communications port settings have to be entered. This is done by pressing the **Setup COM Port** button. Please note that the settings entered here have to match those set in the oscilloscope. The options are:

Port: This is the COM port on the computer that is to be used to communicate with the oscilloscope. The options are COM1 and COM2. since the CD920 controller is normally on port COM1, the oscilloscope is then on port COM2.
Baud rate: This is the data transfer rate, and is normally set to 38400 baud.
Data bits: This is the number of data bits sent per byte transferred and is normally set to 8.
Stop bits: This is the number of stop bits used and is normally set to 1.
Parity: This is the parity convention to be used and is normally set to None.
Flow control: This is the flow control protocol; the value is normally set to Hardware.

The **OK** button is used to accept the new values and to close the form. The **Cancel** button closes the form without accepting the changes.

Setup	×
Settings	
Port	COM2
Baud rate	38400 💌
Data bits	8 💌
Stop bits	1 💌
Parity	None
Flow control	Hardware 💌
	OK Cancel

Figure 3.14: Example of oscilloscope RS-232 settings.

3.5.3 Cryostat

If a cryostat has been attached to the system then pressing the corresponding **Details** button will open the form shown in Figure 3.15. The cryostat uses serial communications to interact with the F900 software.

Cryostat C	onfigura	tion	×
Interface Settin COM Port	gs COM3 💌	Sensor Range Range Lo (K) 10.0	
Baud Rate	9600 💌	Range High (K) 273.0	
GPIB Addr	12 *		
Default		OK Cancel	

Figure 3.15: Example of cryostat settings.

The settings details are:

COM Port: This is the communications port on the computer that is used to talk to the cryostat. The options are COM1, COM2, COM3, COM4. If an oscilloscope has been attached to the system then this setting is normally COM3, otherwise it is normally COM2.

Baud Rate: This is the data transfer rate between the computer and the cryostat and has to be set to the same value on both instruments. The normal setting is 9600 baud.

Range Low: This value is the lower temperature limit of the sensor in Kelvin.

Range High: This value is the upper limit of the sensor in Kelvin.

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

3.5.4 Iris

If a computer-controlled Iris has been installed in the system, then the dialog shown in Figure 3.16 should be opened to set its values. The **High Level Info** grouping contains information on how wide the iris should be

opened. The **Low Level Info** grouping contains factory-set values for the stepper motor controlling the iris. The low-level settings are not normally changed in the field. Details of the settings are:

Description: This value is entered by the user to put a name to the configuration setup. It is only for identification purposes and has no effect on system operation.

Range Low: This is the lowest available opening for the iris (as a percentage).

Range High: This is the highest available opening for the iris (as a percentage.

Initial Position: This is the opening (expressed as a percentage) that the iris will initially move to when the software is started opening for the iris.

Resolution: This is the increment that is added when the iris opening is changed. The control of the iris is such that only integral values can be entered here.

The **Low Level Info** settings should not be altered from their factory settings, without explicit instructions from the Edinburgh Instruments service department. The factory settings are:

Step Multiplier: 1 Step Zero: 0 Step Mode: Half Step Sequence: 2 Ph Step Freq: 4

Iris Co	nfigurat	tion					×
		Description		Default]	
High	Level Info		[Low Level Info			
Rar	ige Low	1		Step Multiplier	1		
Rar	ige High	100		Step Zero	0		
Initia	al Position	1		Step Mode	Ha	lf 💌	
Res	olution	1		Step Sequence	2 F	'n 🔻	
				Step Freq (ms)	4.0	0	
					_		<u> </u>
Defau	lt			OK]	Cancel	

Figure 3.16: Example of iris settings.

3.6 Monochromator Tab

The **Monochromators** tab (shown in Figure 3.17) contains information on the monochromators that have been attached to the system. In a standard configuration there is an Excitation monochromator to select the wavelength of the light entering the sample chamber and an Emission monochromator to select the wavelength of the light exiting the sample chamber. For T geometries there is a second Emission monochromator on the other output arm. For monochromators have more than one grating (three are possible for TM300 and TMS300 monochromators), it is possible to change the selection of grating without going into the details of the selected monochromator.

Detector	s [(CD Controller Li		fetime Cards
Misc.	Lamps	amps Monochromators		Polarisers
Excitation M	ono Use Gr • 1: 2	ating 200.00 - 900.00 nn	n	Details
TM300	Iono Use Gr I: 2	ating 200.00 - 900.00 nn	n	Details
-Emission 2 M	1ono			Details

Figure 3.17: Monochromator tab.

If the monochromator **Details** button is pressed, then the form shown in Figure 3.18 is displayed. This form has several sections, which are:

Present: Whether the monochromator is present on the spectrometer system.

Type: This is the type of the monochromator that is fitted. The options are TM300, TM300, M300, M100. For monochromators with computer-controlled slits (these may be labelled as TMS300 monochromators), select the TM300 and tick the **Present** box in the **Controlled Slits** grouping.

Stepper Details: The details of the stepper motor controlling the monochromator grating(s).

Gratings: The availability of the monochromator gratings is set here and the details for each grating can be displayed.

Wavelength Increments: These are the incremental values that can be used to change the monochromator wavelength in the application software(F900); they are changeable by pressing the **Details** button.

Ports: This grouping is used to set the configuration of the input ports for the Excitation monochromator or the output port for the Emission monochromator(s).

Controlled Slits: If the monochromator has computer-controlled slits fitted, then clicking the **Details** button will allow their settings to be entered.

The settings will now be discussed in greater detail.

Ex Mono Details		×
Present Type: TM300	•	Stepper Details
I: Present 200.00 - 900.00 nm		Details
2: Present No Desc.		Details
🔲 3: Present No Desc.		Details
Wavelength Increments		
0.05,0.10,0.20,0.50,1.00,2.00,5.00		Details
Ports		
Mirror 1 Present	Initial P	osition: Port A 💌
Mirror 2 Present	Initial P	osition: Port A 💌
Controlled Slits		
Present		Details
	OK	Cancel

Figure 3.18: Sample of detailed monochromator settings.

3.6.1 Stepper Details

Em1 Mono MicroStepper Details
Step Period: 4
Step Lower Limit: 0
Step Upper Limit: 8500000
Step Zero: 0
μsteps per full step: 25
Defaults OK Cancel

Figure 3.19: Sample of monochromator micro-stepper settings.

The details for the monochromator stepper motor details are shown in Figure 3.19. These values are normally factory-set and should not be changed without consulting Edinburgh Instruments' service department. The factory-settings are:

Step Period: 4 Step Lower Limit: 0 Step Upper Limit: 8500000 Step Zero: 0 μsteps per full step: 25 A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

3.6.2 Gratings

The **Gratings** grouping allows the user to select what gratings are present. Note that if the monochromator is an M300 or an M100 then only one grating will be available. Pressing the **Details** box for a particular grating opens up the dialog box shown in Figure 3.20.

Ex Grating 1 Details 🛛 🗙
Grating Type: 1800 💌 g/mm
Description: 200.00 - 900.00 nm
Wavelength Data
Range Low: 200.00 🔺 nm Range High: 900.00 🔺 nm
Resolution: 0.05 📥 nm Initial Position: 200.00 🛫 nm
Bandwidth
Linear Dispersion: 1.80 nm/mm Resolution: 0.009 nm
Range Low: 0.02 nm Range High: 27.00 nm
OK Cancel

Figure 3.20: Example of monochromator grating details.

This dialog box allows the user to enter detailed settings for the selected grating as well as display information on the computer-controlled slits attached to the monochromator (if these are fitted). The information that can be entered is:

Grating type: This drop-down box permits the user to enter the groove density (in grooves per mm) of the grating. The allowable values are 50, 75, 150, 300, 600, 840, 1200, 1800, 2400, 3600.
Description: This edit box allows the user to enter a description of the grating. It usually contains the wavelength limits for the grating, although it does not have to. The user should be aware that this description is used in the software to describe the grating, so it is important to enter meaningful text here.

Range Low: This is the lower bandwidth limit of the grating's response function.

Range High: This is the upper bandwidth limit of the grating's response function.

Resolution: This is the limiting resolution of the grating.

Initial Position: This is the position of the grating when the F900 software is initialised.

If the monochromator has computer-controlled slits then the bandwidth limits, resolution, and linear dispersion for the selected grating are shown here. Should the settings for the slits be subsequently changed then these values will also change.

The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

3.6.3 Wavelength Increments

Pressing the **Details** button in the **Wavelength Increments** grouping opens the dialog box shown in Figure 3.21. The step sizes for each step number can be entered, either directly, or by clicking the increment arrows, to change the value of a nominated step. It is normal practice to have the step sizes in order of increasing increments.

MonoSte	pSizes		×
Step Sizes			
	Step 1 (nm)	0.05	
	Step 2 (nm)	0.10	
	Step 3 (nm)	0.20	
	Step 4 (nm)	0.50	
	Step 5 (nm)	1.00	
	Step 6 (nm)	2.00	
	Step 7 (nm)	5.00	
Default		OK Car	icel

Figure 3.21: Example of monochromator step size settings.

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

3.6.4 Ports

If there are mirrors on the entrance (excitation) or exit (emission) monochromators then these should be ticked in the appropriate check box in the **Ports** grouping. Checking Mirror 1 denotes that there is a mirror at the entice port and checking **Mirror 2** indicates that there is a mirror at the output port. One a mirror has been indicated as being present, a drop-down box (**Initial Position**) becomes editable. This box is used to indicate to the F900 software what position the mirror should be in when the software is initialised. The alternatives are Port A and port B.

3.6.5 Controlled Slits

If the monochromator is fitted with computer-controlled slits than this check box should be ticked; otherwise it should be left unticked. Pressing the **Details** button will open the dialog box shown in Figure 3.22. This box allows the user to enter the limiting and initial values for the slits. The details are:

Range Low: This is the smallest slit gap that can be set in the F900 software. The smallest value that can be entered here corresponds to 1 step of the stepper motor, which with a resolution of 200 steps per mm, corresponds to 0.005 mm.

Range High: This is the largest slit gap that can be set in the F900 software. This value will depend on the hardware construction of the slit device; currently this is 15 mm.

Initial Position: This is the slit gap that is set when the F900 software is initialised. **Resolution:** This edit box is used to enter the number of motor steps required to change the slit separation by 1 mm. With existing hardware, this value is 200 steps per mm.

Ex Slit Details	<u> </u>
Slit width data	
Range Low: 0.010 🚔 mm	Initial Position: 0.50 🚔 mm
Range High: 15.00 🚊 mm	Resolution: 200 🚔 steps per mm
Stepper Details	OK Cancel

Figure 3.22: Example of motorised slit settings.

The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box. If the **Stepper Details** button is pressed then the dialog box shown in Figure 3.23 is shown.

🗖 Slit Stepper Details 🛛 🗖 🗙
Step Period: 1.95
Step Zero: 1
Step Multiplier: 1
Step Mode: Full 💌
Step Sequence: 2 Ph 💌
Defaults OK Cancel

Figure 3.23: Example of motorised slit stepper motor details.

This box contains details of the stepper motor settings for the slit device. These values should not normally be changed except by guidance from Edinburgh Instruments service department. The default values for the slit stepper motor are:

Step Period: 1.95 Step Zero: 1 Step Multiplier: 1 Step Mode: Full Step Sequence: 2 Ph

A set of default values can be loaded by pressing the **Default** button. The values are stored by pressing the **OK** button and abandoned by pressing the **Cancel** button. Pressing either of these buttons will close the dialog box.

3.7 Polarisers

The **Polarisers** tab is shown in Figure 3.24. If polarisers have been fitted to the system then the appropriate check boxes should be ticked. The details for a particular polariser can be displayed by pressing the appropriate **Details** button. This will open the dialog box shown in Figure 3.25.

C	D Controller	Life	etime Cards 🛛 🕴
Lamps	Monochroma	tors	Polarisers
			Details
			Details
			Details
		CD Controller Lamps Monochroma	CD Controller Lif Lamps Monochromators

Figure 3.24: Polarisers tab.

The **High Level Info** grouping contains information on the angles that the polariser should be set. The **Low Level Info** grouping contains factory-set values for the stepper motor controlling polariser. The low-level settings are not normally changed in the field. Details of the settings are:

- **Description:** This value is entered by the user to put a name to the configuration setup. It is only for identification purposes and has no effect on system operation.
- **Range Low:** This is the smallest angle that the polariser can be set in the F900 software (expressed in degrees).
- **Range High:** This is the largest angle that the polariser can be set in the F900 software (as a percentage.
- **Initial Position:** This is the initial angle that the polariser will initially move to when the F900 software is initialised.
- **Resolution:** This is the change in value that is added when the polariser angle is incremented in the F900 software..

The **Low Level Info** settings should not be altered from their factory settings, without explicit instructions from the Edinburgh Instruments service department. The factory settings are:

Step Multiplier: not used. Step Zero: 0 Step Mode: Full Step Sequence: Wave Step Freq: 1.60

Excitation Polariser Configuration				
Description Default				
High Level Info		Low Level Info		
Range Low	0.00	Step Multiplier		
Range High	90.00	Step Zero 0		
Initial Position	0.00	Step Mode Full 💌		
Resolution	0.40000	Step Sequence Wave 💌		
		Step Freq (ms) 1.60		
Default		OK Cancel		

Figure 3.25: Example of polariser configuration settings.

4 Software Features

This chapter of the manual describes how to start the F900 software, and how to use the features that are normally visible whenever the software is running. It is assumed that the software has already been installed in the folder 'C:\Program Files\EI\F900'.

With all components of the spectrometer switched on e.g. the Xe900 Xenon lamp and CD920/CD930 controller (switch on the front of the sample chamber); turn on the computer and double-click the F900 icon shown on the desktop. Alternately the F900 program can be reached from going into the 'EI' folder within program files in either windows explorer or in the start menu. Starting the software causes it to detect and initialise all the system components that had been selected in Hardware Configuration.

Important: The software has a protection device (dongle) that needs to be installed in the parallel port LPT1 before starting the software.

4.1 Main Menu and Shortcut Toolbar



Figure 4.1: Main menu and EI toolbar.

The main menu that is displayed at the start up of the F900 software, prior to any data acquisition, is shown in Figure 4.1. This menu comprises drop-down menus for operation of the F900 software and optional shortcut icons for the most commonly used operations. The shortcut toolbar can be enabled/disabled by clicking **View/Tool bar**. The submenu options will be described in the following sections.

The shortcut keys in the toolbar are grouped. The first group of three keys is important for data acquisition. These are the shortcut keys most often used when the user acquires data. The second group of five shortcut keys provide direct access to standard Windows file handling (**Open, Save**), file import and export options and printing facilities. The remainder of the shortcut toolbar are primarily graphics options and data windows options. The functions of the shortcuts are now described.



Print: If a printer is connected directly to your PC, or the PC is networked giving you access to a PC, the button showing a printer symbol enables the user to print a specified measurement container.

Once data has been acquired, it can be viewed in four different ways: 2D, 3D, text, contour. The data plot view buttons enable the user to change the viewing format of data that has been acquired.



A

View As 2D Plot: This option displays the active graph as a two dimensional (2D) plot.



View As 3D Plot: This option displays the active graph as a three dimensional (3D) plot. This option is only available for map-type plots.



View As Text: This option displays the active graph as text.



View As Contour Plot: This option displays the active graph as a contour plot. This option is only available for map-type plots.



Plot options: This button opens a menu that permits the user to change the plot options for the active graph.

The active measurement window can be modified while data is being acquired. The measurement plot can also be altered once the data has been acquired using the following options on the toolbar.

Zoom in: Pressing this button allows the user to zoom into an area of a scan of € particular interest by drawing a rectangle around the chosen area. Zoom out: Pressing this button causes the whole graph to be displayed on screen. Q Cursor on: Pressing this button toggles the visibility of the cursor. The cursor appears 困 as a pair of orthogonal crosshairs, with their intersection being displayed on the plot. Grids on: Pressing this button toggles the visibility of a grid connecting the major tick Ŧ marks on the horizontal and vertical axes. Peaks on: Pressing this button toggles the visibility of the values of the plot maxima and minima. The conditions for setting the peak search criteria can be found in the М Data menu Normalisation: Pressing this button toggles the normalisation the x and y graphs on N the contour plot view. Properties: Clicking this box allows the user to view the instrument and other r settings for a scan. Some of these values may be edited. Join visible: Pressing this button joins all the visible, non-minimised plots into a single graph. This option is only available in 2D view and will only join plots of the same type *****+ as the currently active plot. Split: Pressing this button will separate the plots in the currently active graph into separate graphs, each containing a single plot. Only available in 2D view 2



Extract: Pressing this button will open a dialog box that invites the user to extract a single plot from the currently selected graph.



Logarithmic scale: Pressing this button will display the y-axis as a logarithmic scale and will de-select the **linear scale** button. This option will only work if all the data points have positive y values.



Linear scale: Pressing this button will display the y-axis as a linear scale and will deselect the **logarithmic scale** button.



Abort: Pressing this button stops any measurement mid-scan. A dialog box first appears asking for confirmation of the measurement termination.

4.2 Status Bar

The status bar gives the user an indication of the current mechanical settings of the spectrometer. The status bar can be enabled/disabled by clicking *View/Status Bar*, its layout, for a standard FS920 system, when the software is loaded, is shown in Figure 4.2.

Sam	mple	Excitation 200.00 - 900.00 nm	Emission 1 200.00 - 900.00 nm
1	1	200.00nm, 0.90nm,	200.00nm, 0.90nm,

Figure 4.2: Status Bar of Standard Spectrometer at Start Up

Figure 4.2 shows the wavelength settings of the each monochromator (excitation and emission) and which sample is currently in place (this will always be '1' unless a sample positioner is used, see section 1.2.5).

The information displayed on the status bar alters depending on what the spectrometer is currently undertaking. The bottom left area of the status bar contains information on what process the spectrometer is carrying out at that time. Figure 4.3 shows the status bar as the spectrometer is setting up, before undertaking a spectral measurement. Figure 4.4 shows a status bar while a spectral measurement is being undertaken.

				0 counts
	Repeat	Sample	Excitation	Emission 1 (200.00-900.00nm)
Moving Ex Monochromator to 355.00	1/1	1	500.00nm	650.00nm

Figure 4.3: Status Bar as Spectrometer is Setup for Measurement

				28 counts
Scanning Em1	Repeat	Sample	Excitation	Emission 1 (200.00-900.00nm)
	1/5	1	355.00nm	656.00nm

Figure 4.4: Status Bar During Spectral Measurement

5 Acquiring Spectral Data

5.1 Setup Menu: Signal Rate

The signal rate dialog box (Figure 5.1) is the primary instrumentation setup tool for the F900 software. It allows the signal from the sample under investigation to be optimised before a measurement is performed. The signal rate screen should be used before undertaking any measurements to adjust the following parameters:

- Iris (light levels)
- Wavelength of excitation and emission monochromators.
- Bandwidth of monochromator slits (if computer-controlled slits fitted).
- Polarisation directions (if applicable).
- Excitation source.
- Sample position.
- Detector.

The signal rate screen can also be accessed via the Signal Rate icon on the toolbar.

Signal Rate			? ×
Sample 1			
Excitation Wavelength (n 300.00 🚔 step: 0.50	m) ▼ Δλ: 0.18 ᆍ	Ex Polariser Angle (*): 0.0 -	Source OnF900 © Xe
Em1 Wavelength (nm) 400.00 🚔 step: 0.05	🔹 Δλ: 1.50 🚔	Em1 Polariser	Em1 Detector R955 💌
Iris Setting: 50 🚔			
Temperature: 0.00 °C	🔽 Auto Update	🔽 Display Ref signal	
0 Ref		270 cps	1000
0 Em ⁻		270 cps	1000
		Close	Apply



5.1.1 Dialog Box Settings and Terms

The excitation and emission wavelengths can be set by using the up and down arrows beside the corresponding wavelength edit box. The wavelength increment for each arrow press is determined by the step value, which can beset in the drop-down box adjacent to the wavelength edit box. Alternately, the parameter value can be changed using the keyboard to enter a revised value. Clicking **Apply** or pressing the **Return** key on the keyboard then updates the hardware.

The iris setting can be altered by using the arrows beside the iris setting box. The iris controls the intensity of radiation incident on the sample and is located in the sample chamber just before the reference detector (Figure 1.2). The iris is computer controlled with a stepper motor. The value in the iris box is an integer relating to how much light it lets through. '100' is the maximum setting and corresponds to the iris fully open. '1' is the lowest setting.

If optional polarisers are fitted, the polarisation of the excitation illumination can be set by first moving the excitation polariser to the *"in"* position (into the beam path); clicking the **Ex Polariser** check box and then choosing the required angle in the corresponding drop-down box.

If there is more than one light source connected to the spectrometer, the radio switches beside the names of the sources can be used to toggle between them. Correspondingly, if more than one detector is present, the desired detector can be selected from the drop-down box.

Signal rate dialog boxes show real time readings from the both the reference detector and the emission detector. The reference detector rate can be displayed by ticking the **Display Ref Signal** check box. **Ref** refers to the signal measured on the reference detector; **Em1** refers to the signal measured on the detector connected to the emission monochromator. If a T-Geometry configuration is present, **Em2** refers to the signal measured by the detector connected to the second emission arm. The signal rates are displayed in photon counts per second (cps) and are updated every second. The values are achieved by measuring the detector output for one tenth of a second and multiplying that value by 10. Below the value for the number of counts per second at each detector, there is a rate bar showing the counts with an automatic scaling function. The maximum value of the particular scale is shown to the top right of the rate bar. Although the detectors are used. For analogue detectors (such as an InGaAs detector or the silicon photodiode used as the reference detector), the rate displayed represents a relative irradiative intensity (on an arbitrary scale) measured by the detector.

A temperature probe is fitted to the water cooled sample holder and therefore gives the actual temperature of the sample. The temperature is measured when the signal rate screen is first opened and this measured value is displayed. The **Auto Update** check box facilitates continuous temperature monitoring. When the water cooled sample holder temperature probe is used, the temperature is given in degrees Centigrade (°C), however if an optional cryostat system is fitted, the temperature in the cryostat is displayed in Kelvins (K)

If the optional Sample Positioner is fitted there will be two more tabs present in addition to **Sample 1**. These are called **Sample 2** and **Sample 3**. Each of these tabs contain the same information as **Sample 1** but the settings are 'remembered' so the values for wavelength, polarisation, source, detector and iris can be set up for different samples.

5.1.2 Using the Signal Rate Screen

The main use of the signal rate dialog box is to set up the spectrometer before making a measurement. If the peak excitation and emission wavelengths are known, the monochromators can be set to those wavelengths and then the slits and iris adjusted to set up a bandwidth with sufficient counts to obtain useful spectra. If the peak wavelengths of excitation and emission are not known, the arrows beside the wavelengths can be used to scan through the wavelengths of excitation and emission until a maximum is established. The slits and iris can then be adjusted. Because of an inherent attribute of monochromators with motorised gratings called backlash, for increased accuracy it is best to approach the desired wavelength from the lower wavelength range.

5.2 Scan measurements

The spectral acquisition drop-down menu can be reached through the **New Spectral Measurement** option in the file menu, or by clicking the **Spectral Measurement** icon on the toolbar.

5.2.1 Emission Scan

To perform an emission scan, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down. From this menu select **Emission Scan**. The emission scan dialog box then appears (Figure 5.2). Although it is normal practice to use the **Signal Rate** screen to set up the instrumentation prior to commencing a measurement, some parameters can also be set in the scan setup screens, as will be seen.

Emission Scan Setup ? 🗙
General Excitation Emission1
Position Measurement Mode
Using existing files
Ex Correction Value Measure Simultaneously
Em1 Correction File Data
Emission Scan Parameters
Scan from 360.00 to 450.00 step 10.00 nm
Dwell Time 5.00 Number of Scans 5
Start Cancel Apply

Figure 5.2: Emission scan dialog box.

If the optional sample positioner is being used, the **Position** radio buttons and a group of **Measure** check boxes are available. The **Position** radio buttons allow scans to be set up differently for each sample to be measured. When the measurement is started, the samples measured are dependent on which **Measure** check boxes have been ticked.

If an optional T-Geometry system is being used, **Measurement Mode** radio buttons are present. If **L1** is selected, Emission channel 1 will be used (i.e. data will be acquired from the detector set up to be used on emission channel 1). If **L2** is selected, Emission channel 2 will be used (i.e. data will be acquired from the detector set up to be used on emission channel 2). If **T** is selected, data will be acquired from emission channels 1 and 2 sequentially. Using a **T** measurement mode a wavelength offset can be applied between emission channels 1 and 2. The offset is set on the emission 2 tab (Figure 5.6).

On the General tab select what correction, if any, is to be applied to the data. The correction options are:

Ex Correction Value: This setting corrects for the excitation profile of the spectrometer.
Em1 Correction File Data: This setting corrects for the spectral response of the emission monochromator and detector.
Measure Simultaneously: Checking this setting will mean that the excitation light intensity is measured by the reference detector simultaneously with the emission detector. This option takes account of variations in light source intensity with time as well as spectral variation. Simultaneously measuring the reference detector is quite often a more accurate option, since if the monochromator slit sizes are different from the size they were when the excitation correction file was made, the spectral resolution of the monochromator will alter and therefore the lines of the Xenon spectra measured can be broadened or narrowed.

The standard 450 W Xenon lamp used in the spectrometer is very stable, therefore if correction is required for emission scans, only the **Em1 Correction File Data** check box need be ticked. When any emission correction

is ticked, the software automatically selects the correct file depending on which detector is being used to acquire the data.

On the **Excitation** tab (Figure 5.3), the values for the monochromator wavelength, slit bandwidth, lamp source and polariser that were set in the **Signal Rate** screen are displayed. If desired, the monochromator wavelength, lamp source and optional polarisers, if fitted, can be further modified.

On the **Emission** tab(s) (Figure 5.4), the values for the slit bandwidth, detector and polariser that were set in the **Signal Rate** screen are displayed. If desired, the detector and optional polarisers, if fitted, can be further modified.

In the **Emission Scan Parameters** box (Figure 5.2) enter the wavelength scanning range (**Scan from/to**); step size of scan (*Step*); dwell period at each wavelength (**Dwell Time**) and number of times the scan is to be repeated, and summed (**Number of Scans**).

On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

General Excitation Emission1			
Monochromator Wavelength (nm) 350.00	Polariser In	Angle from	m Vertical (*)
Camp Source	-Bandwidth- Δλ:	0.18	nm



General Excitation Emission1				
	Polariser			
	∏ In	Angle fro	m Vertical (*)	
Detector	Bandwidth			
R955 PCS 1	Δλ:	1.50	nm	

Figure 5.4: Emission scan dialog box, emission tab.
Emission Scan Setup	? X
Position O 1 O 2 O 3	Measurement Mode
Correction Using existing files Ex Correction Value Em1 Correction File Data Em2 Correction File Data	Using reference detector Measure Simultaneously
Emission Scan Parameters	
Scan from 200.00 to 30	00.00 step 1.00 nm
Dwell Time 0.1 Nu	mber of Scans 1
Measure 1 1 2 7 3 Start	Cancel Apply

Figure 5.5: Emission scan dialog box with a T-Geometry and optional sample positioner present.

General Excitation Emission1	Emission2		
Monochromator Em1 to Em2 offset (nm) 20.00	Polariser [] In	Angle fro	om Vertical (°)
Detector InGas PCS Ch3	Bandwidth Δλ:	0.90	nm

Figure 5.6: Emission scan, with emission 2 tab in T-Measurement mode

5.2.2 Excitation Scan

An excitation scan is normally used to obtain excitation spectra of a sample emitting at a specific wavelength. To perform an excitation scan, click the **File** main menu item, then select **New Spectral measurement**. This will display a drop-down. From this menu select **Excitation Scan**. The excitation scan dialog box then appears (Figure 5.7). Although it is normal practice to use the **Signal Rate** screen to set up the instrumentation prior to commencing a measurement, some parameters can also be set in the scan setup screens, as will be seen.

If the optional sample positioner is being used, the **Position** radio buttons and a group of **Measure** check boxes are available. The **Position** radio buttons allow scans to be set up differently for each sample to be measured. When the measurement is started, the samples measured are dependent on which **Measure** check boxes have been ticked.

If an optional T-Geometry system is being used, **Measurement Mode** radio buttons are present. If **L1** is selected, Emission channel 1 will be used (i.e. data will be acquired from the detector set up to be used on

emission channel 1). If **L2** is selected, Emission channel 2 will be used (i.e. data will be acquired from the detector set up to be used on emission channel 2). If **T** is selected, data will be acquired from emission channels 1 and 2 sequentially. Using a **T** measurement mode a wavelength offset can be applied between emission channels 1 and 2. The offset is set on the emission 2 tab.

Excitation Scan Setup ? ×
Position Measurement Mode Image: 1 to 1 t
Correction Using existing files Ex. Correction File Data Em1 Correction Value
Excitation Scan Parameters
Scan from 270.00 to 395.00 step 1.00 nm
Dwell Time 0.20 Number of Scans 5
Measure Start Cancel Apply

Figure 5.7: Excitation scan dialog box.

On the General tab select what correction, if any, is to be applied to the data. The correction options are:

Ex Correction Value: This setting corrects for the excitation profile of the spectrometer. **Em1 Correction File Data:** This setting corrects for the spectral response of the emission monochromator and detector.

Measure Simultaneously: Checking this setting will mean that the excitation light intensity is measured by the reference detector simultaneously with the emission detector. This option takes account of variations in light source intensity with time as well as spectral variation. Simultaneously measuring the reference detector is quite often a more accurate option, since if the monochromator slit sizes are different from the size they were when the excitation correction file was made, the spectral resolution of the monochromator will alter and therefore the lines of the Xenon spectra measured can be broadened or narrowed.

The standard 450 W Xenon lamp used in the spectrometer is very stable, therefore if correction is required for emission scans, only the **Em1 Correction File Data** check box need be ticked. When any emission correction is ticked, the software automatically selects the correct file depending on which detector is being used to acquire the data.

On the **Excitation** tab (Figure 5.8), the values for the monochromator wavelength, slit bandwidth, lamp source and polariser that were set in the **Signal Rate** screen are displayed. If desired, the monochromator wavelength, lamp source and optional polarisers, if fitted, can be further modified.

General Excitation Emission1		
	Polariser	Angle from Vertical (*)
	I IN	
Lamp Source	Bandwidth	
● Xe900 ○ uFlamp	Δλ:	5.00 nm

Figure 5.8: Excitation scan dialog box, excitation tab.

On the **Emission** tab(s) (Figure 5.9), the values for the slit bandwidth, detector and polariser that were set in the **Signal Rate** screen are displayed. If desired, the detector and optional polarisers, if fitted, can be further modified.

General Excitation Emission1		
Monochromator Wavelength (nm) 592.30	Polariser	Angle from Vertical (*)
Detector R955 PCS 1	Bandwidth Δλ:	1.50 nm

Figure 5.9: Excitation scan dialog box, emission tab.

In the **Excitation Scan Parameters** box (Figure 5.7) enter the wavelength scanning range (**Scan from/to**); step size of scan (*Step*); dwell period at each wavelength (**Dwell Time**) and number of times the scan is to be repeated, and summed (**Number of Scans**).

On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

5.2.3 Synchronous Scan

In synchronous scans, both excitation and emission monochromators are scanned synchronously with a preset offset. These scans can be used in order to separate and identify fluorescence components in mixtures. For low concentrations, each fraction of the molecules has a peak at a characteristic wavelength of the synchronous spectrum.

To perform a synchronous scan, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down. From this menu select **Synchronous Scan**. The excitation scan dialog box then appears (Figure 5.10). Although it is normal practice to use the **Signal Rate** screen to set up the instrumentation prior to commencing a measurement, some parameters can also be set in the scan setup screens, as will be seen.

Synchronous Scan Setup 🛛 🔹 🗙
General Excitation Emission1
Position Measurement Mode © 1 C 2 C 3
Correction Using existing files Ex. Correction File Value Em1 Correction File Data Using reference detector Measure Simultaneously
Excitation Scan Parameters
Scan from 300.00 to 400.00 step 1.00 nm
Dwell Time 0.20 Number of Scans 1
Measure Start Cancel Apply

Figure 5.10: Synchronous scan dialog box.

If the optional sample positioner is being used, the **Position** radio buttons and a group of **Measure** check boxes are available. The **Position** radio buttons allow scans to be set up differently for each sample to be measured. When the measurement is started, the samples measured are dependent on which **Measure** check boxes have been ticked.

If an optional T-Geometry system is being used, **Measurement Mode** radio buttons are present. If **L1** is selected, Emission channel 1 will be used (i.e. data will be acquired from the detector set up to be used on emission channel 1). If **L2** is selected, Emission channel 2 will be used (i.e. data will be acquired from the detector set up to be used on emission channel 2). If **T** is selected, data will be acquired from emission channels 1 and 2 sequentially. Using a **T** measurement mode means that synchronous scans with two different offset wavelengths can be acquired at the same time.

On the General tab select what correction, if any, is to be applied to the data. The correction options are:

Ex Correction Value: This setting corrects for the excitation profile of the spectrometer. **Em1 Correction File Data:** This setting corrects for the spectral response of the emission monochromator and detector.

Measure Simultaneously: Checking this setting will mean that the excitation light intensity is measured by the reference detector simultaneously with the emission detector. This option takes account of variations in light source intensity with time as well as spectral variation. Simultaneously measuring the reference detector is quite often a more accurate option, since if the monochromator slit sizes are different from the size they were when the excitation correction file was made, the spectral resolution of the monochromator will alter and therefore the lines of the Xenon spectra measured can be broadened or narrowed.

The standard 450 W Xenon lamp used in the spectrometer is very stable, therefore if correction is required for emission scans, only the **Em1 Correction File Data** check box need be ticked. When any emission correction is ticked, the software automatically selects the correct file depending on which detector is being used to acquire the data.

On clicking Start, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

On the **Excitation** tab (Figure 5.11), the values for the monochromator wavelength, slit bandwidth, lamp source and polariser that were set in the **Signal Rate** screen are displayed. If desired, the monochromator wavelength, lamp source and optional polarisers, if fitted, can be further modified.

General Excitation Emission1			
	Polariser		V / 100
] In	Angle from	m Vertical (*)
Lamp Source	Bandwidth		
C uFlamp	Δλ:	5.00	nm

Figure 5.11: Synchronous scan excitation tab.

On the **Emission** tab(s)) (Figure 5.12), the values for the slit bandwidth, detector and polariser that were set in the **Signal Rate** screen are displayed. If desired, the detector and optional polarisers, if fitted, can be further modified.

General Excitation Emission1		
Monochromator	Polariser	
Ex to Em1 offset (nm)	🗖 In	Angle from Vertical (*)
10.00		0.0 💌
Detector	Bandwidth	
R928-P	Δλ:	5.00 nm
C R5509		

Figure 5.12: Synchronous scan emission tab.

In the Excitation Scan Parameters box (Figure 5.10) enter the wavelength scanning range (Scan from/to); step size of scan (*Step*); dwell period at each wavelength (Dwell Time) and number of times the scan is to be repeated, and summed (Number of Scans).

On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

5.2.4 Kinetic Scan

A kinetic scan monitors the luminescence intensity as a function of time subsequent to a single pulse of light from the Xe900 lamp. The lamp is shuttered on for a user-defined period of time and the decay in the luminescent intensity, subsequent to this excitation pulse, is monitored over a user-defined period of time.

To perform a kinetic scan, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down. From this menu select **Kinetic Scan**. The excitation scan dialog box then appears (Figure 5.13).

If the optional sample positioner is being used, the **Sample** radio buttons and a group of **Measure Samples** check boxes are available. The **Sample** radio buttons allow scans to be set up differently for each sample to be measured. When the measurement is started, the samples measured are dependent on which **Measure Samples** check boxes have been ticked.

Kinetic Scan Setup	? ×
Sample © 1 © 2 © 3	Measure Samples
Sample Set Up	
Scan Time (s) 300.0	No. of Data Points: 301
Use Shutter	seconds for 8.0 🔹 seconds
Start	Cancel Apply

Figure 5.13: Kinetic scan setup dialog box.

The settings that can be made and viewed are:

Scan Time: This is the duration of a scan. The value can be entered directly into the edit box or adjusted using the arrows at the side of the box. The time increment for each arrow click is the value of the **Time Resolution** edit box. Changes made with the arrows are implemented immediately, however directly-entered txt is only implemented when the **Apply** button has been pressed.

Time Resolution: This is the time interval between consecutive data acquisitions. This value can be entered directly into the edit box or adjusted using the arrowed buttons. The time increment for each arrow click is 0.1 s. Changes made with the arrows are implemented immediately, however directly-entered text is only implemented when the **Apply** button has been pressed.

No. of Data Points: The number of data points is calculated by dividing the **ScanTime** by the **Time Resolution**. It is updated immediately if changes have been made with arrow buttons or, alternatively, after the **Apply** button has been pressed.

Use Shutter: Checking this box will enable the shutter in front of the Xe lamp.

Open shutter after: This is the delay after the start of the experiment that the shutter in front of the Xe lamp is opened. This value can be entered directly into the edit box or adjusted using the arrowed buttons. The time increment for each arrow click is 1.0 s. Changes made with the arrows are implemented immediately, however directly-entered text is only implemented when the **Apply** button has been pressed.

for: This is the time that the shutter will be open. As before, this value can be entered directly into the edit box or adjusted using the arrowed buttons. The time increment for each arrow click is 1.0 s. Changes made with the arrows are implemented immediately, however directly-entered text is only implemented when the **Apply** button has been pressed.

Any directly-entered text in the text boxes will be implemented when the **Apply** button is pressed. On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied. The dialog box is closed by pressing the **Cancel** button.

Sample Set Up: Pressing this button will permit the setting of the parameters for the sample currently selected in the **Sample** radio group. The dialogs for this purpose are shown in Figure 5.16.

ample Set up 🛛 🔋 🗙
Set up Sample © 1
Mode Excitation Emission 1
Measurement Mode
OK Cancel
(a)

Mode Excitation E	mission 1	
Wavelength (nm) 300.00 - Polariser Angle: 0.0 -	Increment (nm) 0.05 • Light Source • nF900 • Xe	
Iris Setting: 5 🛬	Δλ: 0.18	nm
(b))	
Mode Excitation E	mission 1	
Wavelength (nm) 400.00 + Polariser Angle: 0.0 +	Increment (nm) 0.05 • Detector • R955 • PCS 1	
	Δλ: 1.50	nm

Figure 5.14: Kinetic scan sample set up dialogs with (a) Mode tab, (b) Excitation tab and (c) Emission tab shown.

There are three instrumentation tabs, which are:

Mode: This tab is shown in Figure 5.14(a) and it permits the user to select the measurement mode, based on the system geometry, for the measurement.

Excitation: This tab is shown in Figure 5.14(b) and it permits the user to modify the instrumentation associated with the excitation monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The Light Source radio group is used to select the source to be used in the experiment.
- d. If a computer-controlled iris is fitted, the opening of the iris (as a percentage) can be set using the **Iris Setting** dialog box.

e. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

Emission: This tab is shown in Figure 5.14(c) and it permits the user to modify the instrumentation associated with the emission monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The Detector radio group is used to select the detector to be used in the experiment.
- d. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

5.2.5 Cryostat Temperature Scan

This measurement requires an Oxford cryostat and temperature controller and these settings need to be made in the hardware configuration (Chapter 3) before the menu item for a temperature scan become available. The instructions for set-up and use of a cryostat may be found in the cryostat manuals supplied. A temperature scan measures the fluorescence of a sample at fixed excitation and emission wavelengths over a temperature range.

To perform a temperature scan, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down. From this menu select **Temperature Scan**. The temperature scan dialog box then appears (Figure 5.15). Although it is normal practice to use the **Signal Rate** screen to set up the instrumentation prior to commencing a measurement, some parameters can also be set in the scan setup screens, as will be seen.

During the temperature scan, the cryostat takes the temperature to the lowest point in the temperature range. Once the cryostat has reached the temperature, it has to stay within a preselected temperature range for a given time to make sure of its stability, the emission is then measured for a given time. The temperature in the cryostat sample holder then increases to the next step in the scan. These settings are altered in the **Temperature Scan Parameters** box (Figure 5.15). The parameters that can be entered here are:

Scan From/To/Step: Enter the temperature range of interest in Kelvin (Scan from/to) and the number of degrees Kelvin between each measurement (Step). Stabilisation Time: Enter the time that the cryostat has to stay at each temperature before a measurement is acquired.

Accuracy Band: Enter the temperature tolerance range (in Kelvin) within which is the temperature is assumed to be stable.

Dwell Time: Enter the time data is acquired at each wavelength in seconds.

Auto PID: The Proportional, Integral and Derivative settings can either be set manually on the cryostat controller or selected automatically from the learnt characteristics of the system by means of the Auto PID check box in the **Temperature Scan Parameters**. When the cryostat temperature is set to go to a higher temperature, the temperature increase has a hysteresis profile. This is one of the reasons for a stabilisation time. In order to minimise the hysteresis while maximising the speed of the temperature increase, the proportional, integral and derivative settings can be altered. We recommend that the temperature controller automatically sets these. More information on this can be found in the cryostat instruction manuals.

Temperature Scan Setup (Cryostat) 💦 🛛 🔋 🗙
General Excitation Emission1
Position Measurement Mode
Using existing files
Ex Correction File Value 🔲 Measure Simultaneously
Em1 Correction File Value
Scan from 100.00 to 150.00 step 1.00 K
Accuracy Band ± 0.10 Stabilisation Time 1
Dwell Time 0.10
Auto PID
Start Cancel Apply

Figure 5.15: Temperature Scan Setup Dialog box

If the optional sample positioner is being used, the **Position** radio buttons and a group of **Measure** check boxes are available. The **Position** radio buttons allow scans to be set up differently for each sample to be measured. When the measurement is started, the samples measured are dependent on which **Measure** check boxes have been ticked.

If an optional T-Geometry system is being used, **Measurement Mode** radio buttons are present. If **L1** is selected, Emission channel 1 will be used (i.e. data will be acquired from the detector set up to be used on emission channel 1). If **L2** is selected, Emission channel 2 will be used (i.e. data will be acquired from the detector set up to be used on emission channel 2). If **T** is selected, data will be acquired from emission channels 1 and 2 sequentially. Using a **T** measurement mode a wavelength offset can be applied between emission channels 1 and 2. The offset is set on the emission 2 tab.

On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

On the **General** tab, select the required measurement mode and the correction to be applied to the data. Since the excitation and emission wavelengths are fixed, the spectral correction files need not be used. To correct for any variation in lamp intensity over the course of the measurement, the **Measure Simultaneously** box should be checked. Checking this setting will mean that the excitation light intensity is measured by the reference detector simultaneously with the emission detector. This option takes account of variations in light source intensity with time as well as spectral variation.

General Excitation Emission1	1		
Monochromator	Polariser		
Wavelength (nm)	🗖 In	Angle fro	om Vertical (*)
200.00		0.00	-
Lamp Source	Bandwidth		
Xe900	Δλ:	5.00	nm
C uFlamp			

Figure 5.16: Temperature Scan Excitation Tab

On the **Excitation** tab (Figure 5.16), the values for the monochromator wavelength, slit bandwidth, lamp source and polariser that were set in the **Signal Rate** screen are displayed. If desired, the monochromator wavelength, lamp source and optional polarisers, if fitted, can be further modified.

General Excitation Emission1			
Monochromator	Polariser	Angle from) (ortical (°)
200.00	1_] m		
Detector	-Bandwidth-		
R928-P R5509	Δλ:	5.00	nm
1 1 1 2 2 2 2			

Figure 5.17: Temperature Scan Emission Tab

On the **Emission** tab(s) (Figure 5.17), the values for the slit bandwidth, detector and polariser that were set in the **Signal Rate** screen are displayed. If desired, the detector and optional polarisers, if fitted, can be further modified.

5.3 Correction Scans

Spectral correction is necessary to obtain the true excitation and emission spectra of the sample, free from any instrumental effects. Comprehensive spectral correction, using previously measured correction files, is standard practice when using fluorescence spectrometers.

5.3.1 Emission Correction Scan

Raw emission spectra are affected by the monochromator efficiency and the spectral response of the detector. Unique correction files for each spectrometer are obtained during calibration at the Edinburgh Instruments factory using calibrated light sources. However, if the need arises to measure additional emission correction files (for instance, if a new detector is installed onto the system) you must firstly obtain the appropriate calibrated sources from Edinburgh Instruments Ltd. Once you are in receipt of these, proceed in the following manner to measure a new emission correction file.

Attach the calibration source to the side of the sample chamber at 90° to the emission channel. Place a highly scattering object (such as a PTFE scatterer) in the sample holder at 45° to the calibration lamp and emission monochromator.

Emission1 Correctio	n Scan Setup	? X
Emission1 Lamp Source	Polariser	
tungsten	☐ [n] Angle from Ve 0.0	rtical (*)
Detector	Bandwidth	
R928-P R5509	Δλ: 0.90 nn	n
Emission Scan Parameters		
Scan from 600.00	to 900.00 step 1.00	nm
Dwell Time 0.20	Number of Scans 5	
	Start Cancel	Apply

Figure 5.18: Emission Correction Dialog Box

To perform an emission correction scan, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down menu. From this menu, select **Correction Scan** and on the popup menu click the required emission arm to be measured. The emission correction dialog box will be opened (Figure 5.18). The following items can be edited:

Lamp Source: Enter a description for the calibration source (e.g. Tungsten or Deuterium).

Polariser: If polarisation is to be used then check the **In** box and select the **Angle from Vertical** from the available options.

Detector: Select a detector from the options available.

Scan from/to/step: Enter the wavelength scanning range and the size of the step.

Dwell Time: Enter the dwell period at each wavelength.

Number of Scans: Enter the number of times that the scan has to be repeated and summed.

On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

As the calibration source has its own spectral profile, the data measured has to be divided by this profile before it can be used as a correction file.

Tip:

The spectral range of an emission correction file may be extended by measuring the response of two different calibration sources (e.g. Tungsten and Deuterium), dividing the measurement with their source profiles and then appending the resulting sets of data.

5.3.2 Excitation Correction Scans

Uncorrected excitation spectra are effected by the spectral output of the light source and the throughput of the monochromator. The correction file is obtained by using the built-in calibrated reference detector that monitors a fraction of the excitation light. This scan requires specific settings to be made for the monochromator. To set these, go to the **Signal Rate** menu and enter the following values:

- a) Set the excitation wavelength to 473 nm.
- b) Set the spectral resolution of the slits to 1 nm (equivalent to 0.55 mm for a single monochromator with an 1800 grooves/mm grating).

Excitation (Ref.) Cor	rection	Scan Se	etup ? 🗙
Excitation (Her.)	Polariser		
	🗆 ln	Angle from	m Vertical (*)
		0.0	•
Lamp Source	Bandwidth		
• Xe900	Δλ:	0.90	nm
Excitation Scan Parameters			
Scan from 200.00 t	o 900.00	step 1.0	0 nm
Dwell Time 0.20	Number of	Scans 5	
	Start	Cancel	Apply

Figure 5.19: Excitation Correction Dialog Box

To perform an emission correction scan, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down menu. From this menu, select **Correction Scan** and on the popup menu click **Ex (Ref. Detector)** The excitation correction scan setup dialog box is then opened (Figure 5.19). The following items can be edited:

Lamp Source: Select the required excitation source (normally Xe Lamp).

Polariser: If required, create correction scans for various excitation polarisations can be created, but for the standard correction file, the polariser is left out of the beam path.

Scan from/to/step: Enter the wavelength scanning range and the size of the step. These values should normally be: **From** = 200 nm, **To** = 900 nm, **Step** = 1 nm (for a 1800 grooves/mm grating blazed at 500 nm).

Dwell Time: Enter the dwell period at each wavelength. This should normally be set to 0.2 s.

Number of Scans: Enter the number of times that the scan has to be repeated and summed. This should normally be set to 5.

On clicking Start, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

Once the scan has been completed it should be saved in folder location C:\Program Files\EAI\Correction.

In order to install the correction file so that it is used by the software, follow these steps:

- a. On the F900 main menu bar go to Options/Correction Files.
- b. Go to the Excitation tab and click the Modify button.
- c. Using the open dialog box, locate the correction file you saved, i.e. in C:\Program Files\EAI\Correction.
- d. The path to the correction file should now appear below 'Ex Correction File', if so click Ok
- e. The software should now be configured to work with the new correction file.

<u>Notes</u>

The Silicon reference detector should be limited to 200 -1100 nm.

The excitation correction file should be repeated periodically and when a bulb is changed, as pressure broadening can occur with the Xenon spectral lines.

5.4 Map measurements

5.4.1 Emission Map

Map measurements consist of a pre-defined set of basic scans, each done at a different physical attribute. An emission map performs a set of emission scans for a set of pre-defined excitation wavelengths.

Emission Map Setup	? ×
General Excitation Emission1	,
Position	Measurement Mode
• 1 • C 2 • C 3	© L1 O L2
Correction Using existing files	Using reference detector
Ex Lorrection Value Em1 Correction File Data	
Emission Map Parameters	
Ex scan from 270.00 to 3	395.00 step 10.00 nm
Em scan from 365.00 to 5	500.00 step 10.00 nm
Dwell Time 0.10 N	umber of Scans 1
Measure	t Cancel Apply

Figure 5.20: Emission Map Dialog Box with general tab shown.

To perform an emission map, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down. From this menu select **Emission Map**. The emission map dialog box then appears (Figure 5.20). Although it is normal practice to use the **Signal Rate** screen to set up the instrumentation prior to commencing a measurement, some parameters can also be set in the scan setup screens, as will be seen.

If the optional sample positioner is being used, the **Position** radio buttons and a group of **Measure** check boxes are available. The **Position** radio buttons allow scans to be set up differently for each sample to be

measured. When the measurement is started, the samples measured are dependent on which **Measure** check boxes have been ticked.

General Excitation Emission1		
	Polariser	Angle from Vertical (*)
Camp Source	Bandwidth Δλ:	5.00 nm

Figure 5.21: Excitation tab for an emission map.

General Excitation Emission1			
I Fin1 start > Ex Start offset (nm)	Polariser	Angle fror	n Vertical (*)
10.00		0.0	•
Detector	Bandwidth		
R928-P R5509	Δλ:	5.00	nm

Figure 5.22: Emission tab for an emission map.

On the General tab select what correction, if any, is to be applied to the data. The correction options are:

Ex Correction Value: This setting corrects for the excitation profile of the spectrometer. **Em1 Correction File Data:** This setting corrects for the spectral response of the emission monochromator and detector.

Measure Simultaneously: Checking this setting will mean that the excitation light intensity is measured by the reference detector simultaneously with the emission detector. This option takes account of variations in light source intensity with time as well as spectral variation. Simultaneously measuring the reference detector is quite often a more accurate option, since if the monochromator slit sizes are different from the size they were when the excitation correction file was made, the spectral resolution of the monochromator will alter and therefore the lines of the Xenon spectra measured can be broadened or narrowed.

The standard 450 W Xenon lamp used in the spectrometer is very stable, therefore if correction is required for emission scans, only the **Em1 Correction File Data** check box need be ticked. When any emission correction is ticked, the software automatically selects the correct file depending on which detector is being used to acquire the data.

In the Emission Map Parameters box, the following parameters can be entered:

Ex scan from/to/step: Enter the excitation wavelength scanning range and step size.

Em scan from/to/step: Enter the emission wavelength scanning range and step size.

Dwell Time: Enter the dwell time at each wavelength.

Number of scans: Enter the number of times that the scan is to be repeated and summed.

The number of emission scans that will be completed in order to construct the map will be the number of steps in the excitation wavelength range.

On the **Excitation** tab (Figure 5.21), the values for the slit bandwidth, lamp source and polariser that were set in the **Signal Rate** screen are displayed. If desired, the lamp source and optional polarisers, if fitted, can be further modified.

On the **Emission** tab(s) (Figure 5.22), the values for the slit bandwidth, detector and polariser that were set in the **Signal Rate** screen are displayed. If desired, the detector and optional polarisers, if fitted, can be further modified. By default, if the scan ranges of the excitation and emission monochromators overlap then it is possible for the excitation wavelength to be greater than the emission wavelength. If this is not desired then the **Em1 start > Ex** box should be checked and a value entered in the **Start offset** edit box. The effect of this will be to override the settings in the **Emission Map Parameters** box by ensuring that the emission wavelength starts at the offset value greater than the excitation wavelength.

On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

5.4.2 Synchronous Map

To perform a synchronous map, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down. From this menu select **Synchronous Map**. The synchronous map dialog box then appears (Figure 5.23). Although it is normal practice to use the **Signal Rate** screen to set up the instrumentation prior to commencing a measurement, some parameters can also be set in the scan setup screens, as will be seen.

If the optional sample positioner is being used, the **Position** radio buttons and a group of **Measure** check boxes are available. The **Position** radio buttons allow scans to be set up differently for each sample to be measured. When the measurement is started, the samples measured are dependent on which **Measure** check boxes have been ticked.

On the General tab select what correction, if any, is to be applied to the data. The correction options are:

Ex Correction Value: This setting corrects for the excitation profile of the spectrometer. **Em1 Correction File Data:** This setting corrects for the spectral response of the emission monochromator and detector.

Measure Simultaneously: Checking this setting will mean that the excitation light intensity is measured by the reference detector simultaneously with the emission detector. This option takes account of variations in light source intensity with time as well as spectral variation. Simultaneously measuring the reference detector is quite often a more accurate option, since if the monochromator slit sizes are different from the size they were when the excitation correction file was made, the spectral resolution of the monochromator will alter and therefore the lines of the Xenon spectra measured can be broadened or narrowed.

The standard 450 W Xenon lamp used in the spectrometer is very stable, therefore if correction is required for emission scans, only the **Em1 Correction File Data** check box need be ticked. When any emission correction is ticked, the software automatically selects the correct file depending on which detector is being used to acquire the data.

Synchronous Map Setup ? 🗙
General Excitation Emission1
Position • 1 • 2 • 3 • 1 • L2
Correction Using existing files Ex Correction File Value Em1 Correction File Data Using reference detector Measure Simultaneously
Synchronous Map Parameters
Ex scan from 270.00 to 395.00 step 1.00 nm
Em offset from 0.00 to 500.00 step 1.00 nm
Dwell Time 0.10 Number of Scans 1
Measure Measure Cancel Apply

Figure 5.23: Synchronous map dialog box with general tab shown.

In the Emission Map Parameters box, the following parameters can be entered:

Ex scan from/to/step: Enter the excitation wavelength scanning range and step size.

Em scan from/to/step: Enter the emission wavelength scanning range and step size.

Dwell Time: Enter the dwell time at each wavelength.

Number of scans: Enter the number of times that the scan is to be repeated and summed.

The number of emission scans that will be completed in order to construct the map will be the number of steps in the excitation wavelength range.

On the **Excitation** tab (Figure 5.24), the values for the slit bandwidth, lamp source and polariser that were set in the **Signal Rate** screen are displayed. If desired, the lamp source and optional polarisers, if fitted, can be further modified.

On the **Emission** tab(s) (Figure 5.25), the values for the slit bandwidth, detector and polariser that were set in the **Signal Rate** screen are displayed. If desired, the detector and optional polarisers, if fitted, can be further modified.

On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

General Excitation Emission1		
	Polariser	
	🗖 In	Angle from Vertical (*)
		0.0 💌
Lamp Source	Bandwidth	
	Δλ:	5.00 nm
C uFlamp		



General Excitation Emission1			
	Polariser		
	🗖 In	Angle fro	m Vertical (*)
		0.0	•
Detector	Bandwidth		
R928-P	Δλ:	5.00	nm
C R5509			

Figure 5.25: Emission tab for a synchronous map.

5.4.3 Temperature Map

The temperature map scan requires an Oxford cryostat and temperature controller to be present. To perform a temperature map, click the **File** main menu item, then select **New Spectral measurement**. This will display a drop-down. From this menu select **Temperature Map**. The temperature map dialog box then appears (Figure 5.26).

The parameters that can be set are:

Scan Type: Select the type of scan to be performed from the selection of radio butoons available. The scan can then be set up by pressing the ... **Scan Setup** button. This will open the menus described in Section 5.2.1 (for emission scans), Section 5.2.2 (for excitation scans) or Section 5.2.3 (for synchronous scans).

Scan From/To/Step: Enter the temperature range of interest in Kelvin (Scan from/to) and the number of degrees Kelvin between each measurement (Step).

Stabilisation Time: Enter the time that the cryostat has to stay at each temperature before a measurement is acquired.

Accuracy Band: Enter the temperature tolerance range (in Kelvin) within which is the temperature is assumed to be stable.

Auto PID: The Proportional, Integral and Derivative settings can either be set manually on the cryostat controller or selected automatically from the learnt characteristics of the system by means of the Auto PID check box in the **Temperature Scan Parameters**. When the cryostat temperature is set to go to a higher temperature, the temperature increase has a hysteresis profile. This is one of the reasons for a stabilisation time. In order to minimise the hysteresis while maximising the speed of the temperature increase, the

proportional, integral and derivative settings can be altered. We recommend that the temperature controller automatically sets these. More information on this can be found in the cryostat instruction manuals.

Temperature Map Setup 🛛 🛛 🛛 🖓 🗙	
Scan Type © Emission Scane © Synchronous Scan	
Emission Scan Setup	
Temperature Scan Parameters	
Scan from 295.0 to 297.0 step 1.0 K	
Accuracy Band ± 0.2 Stabilisation Time 5	
Auto PID	
Start Cancel Appl	у

Figure 5.26: Temperature map dialog box.

On clicking **Start**, all sample values are rounded to their hardware device resolution, then checked against their legal range. Any illegal values will be notified and a scan cannot start until these have been remedied.

Select the required cryostat control mode when prompted.

5.5 Anisotropy measurements (performed using optional polarisers)

If optional, computer controlled polarisers (P920) are present, anisotropy scans for both emission scans and excitation scans can be performed. When fluorescent samples are excited by polarised light, photoselection of the fluorophores results, depending on their orientation relative to the direction of polarisation of the light.

Principally, an anisotropy calculation can be made without G-factor correction by loading the .VV and .VH files only. However, because the G-factor is dependent on the emission wavelength, the user should apply the G-factor correction to the fluorescence anisotropy calculation. The G-factor (which is a G-factor file in steady state spectroscopy) is calculated from the .HV and .HH files as follows:

$$G_i = \frac{I_{HHi}}{I_{HVi}}$$

The steady state fluorescence anisotropy is then calculated as follows:

$$\mathbf{r_i} = \frac{\mathbf{G_i} \cdot \mathbf{I_{VVi}} - \mathbf{I_{VHi}}}{\mathbf{G_i} \cdot \mathbf{I_{VVi}} + 2 \cdot \mathbf{I_{VHi}}}$$

5.5.1 Emission Anisotropy Scan

To perform an emission scan, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down. From this menu select **Anisotropy Scan**. From the pull-down menu select **Emission**. The emission anisotropy scan dialog box then An emission anisotropy scan is set up in the same manner as a normal emission scan. Please see Section 5.2.1 for instructions on how to do this. Additionally there is a tab containing settings for the anisotropy measurements (Figure 5.27).

General Excitation Emission1	Anisotropy
Ex Polariser I⊄ Ex Vertical	Em1-Em2 Polarisation C Vertical Horizontal C V then H Angle (*) O V then H V then H
 G Factor Measure curves HH and H Prompt for previously measure 	V and ratio ured G Factor curve on completion

Figure 5.27: Anisotropy tab for an emission scan.

The anisotropy dialogue box allows the user to set up the procedure used to make an anisotropy measurement in the measurement screen [Measure - Anisotropy].

For measurements in **L** geometry only the two excitation polariser options and one emission polariser option (V then H) is active.

If the vertical polarisation direction is selected for the excitation channel the spectrally resolved fluorescence intensities I_{VV} and I_{VH} will be sequentially measured. If the horizontal polarisation direction has been selected the excitation scans I_{HV} and I_{HH} , required for G-factor correction, will be sequentially measured

In **T-geometry**, six different modes of automatic polarised fluorescence measurements can be set up, with three of them used to measure curves in order to determine the G-factor.

The anisotropy set-up dialogue box sets up anisotropy measurements to be measured in the measurement screen **[Measure - Anisotropy]**.

If the vertical polarisation direction is selected for the excitation channel and the first of the three options is selected for the emission channels then I_{VV} and I_{VH} will be respectively and simultaneously measured at the two emission arms.

If the vertical polarisation direction is selected for the excitation channel and the second of the 3 options is selected for the emission channels I_{VV} and I_{VH} will be measured at emission arm 1 subsequently, whereas a user defined polarised measurement will be made simultaneously in the second emission arm. If the vertical polarisation direction is selected for the excitation channel and the last of the 3 options is selected for the emission channels the spectrally resolved fluorescence intensities I_{VV} will be measured at both emission arms simultaneously and I_{VH} will be measured for both arms afterwards.

The measurements for horizontal oriented polariser in the excitation channel are similar to the measurements described above. They are used to make the G factor correction.

All polarised measurements are stored automatically with the file extensions ".VV", ".VH", ".HV", and ".HH", dependent upon the orientation of the excitation and emission polarisers.

5.5.2 Excitation Anisotropy Scan

To perform an excitation anisotropy scan, click the **File** main menu item, then select **New Spectral measurement.** This will display a drop-down. From this menu select **Anisotropy Scan**. From the pull-down menu select **Excitation**. The excitation anisotropy scan dialog box then appears An excitation anisotropy scan is set up in the same manner as a normal emission scan. Please see Section 5.2.2 for instructions on how to do this. Additionally there is a tab containing settings for the anisotropy measurements (Figure 5.28).

General Excitation Emission1	Anisotropy
Ex Polariser	Em1-Em2 Polarisation C Vertical Horizontal C V then H Angle (*) O V then H V then H
 ✓ G Factor ○ Measure at peak ⊙ Use value 1.0000 	

Figure 5.28: Anisotropy tab for an excitation scan.

For an excitation scan (where the G-factor file should be constant) the G-factor can be modified in the **G Factor** box.

6 Acquiring Lifetime Data

If the spectrometer system is fitted with a TCC900 card the lifetime measurements can be performed. As with spectral measurements, the **Signal Rate** dialog box is used to prepare the instrumentation setups for an experiment. Details of the **Signal Rate** settings can be found in Section 5.1. Once the hardware settings have been completed, the measurement parameters can be set by selecting the **New Lifetime Measurement** menu item in the **File** header, or by pressing the **Lifetime Measurement** toolbar button. This will open the following options:



Figure 6.1: Lifetime measurement options

These options will now be described in greater detail.

6.1 Manual Lifetime

When Manual Lifetime is selected then a dialog box similar to the one shown in Figure 6.2 will open:

Manual	measur	ement	? ×
▶ <u>N</u> ew	∎ <u>P</u> ause	■ <u>S</u> top	Window New
⊂SetupSam ●1 C	ple 2 O 3	🗖 IRF	O Add
Mode E	xcitation E	mission 1	1
-Measure	.1 O I	L2 O	Г
🗌 Live	Clos	:e	Apply

Figure 6.2: Manual lifetime dialog box with *Mode* geometry options selected.

This box permits the user to:

a. Start a measurement by pressing the New button.

?

Window

New

O Add

nm

×

- b. Temporarily halt a measurement sequence by pressing the Pause button.
- c. Permanently halt a measurement by pressing the Stop button.
- d. Choose whether data should be displayed in the currently active data window (**Add**) or be shown in a new window (**New**).
- e. For systems with multiple sample chamber positions, the desired position can be selected using the **Set up Sample** radio group.
- f. If the purpose of the measurement is to obtain the system response, without any sample present, then the **IRF** checkbox can be ticked to indicate that the data can be used in deconvolution to remove the effect of the measurement system from acquired sample data.
- g. If the **Live** check box is ticked then monochromator wavelengths can be changed during the course of a measurement, otherwise they are only changed at the beginning of a measurement.

Pressing the **Apply** button makes any changes effective and pressing the **Close** button will close the dialog box.



Figure 6.3: Manual lifetime dialog box with *Excitation* monochromator options selected.

Figure 6.4: Manual lifetime dialog box with *Emission* monochromator options selected.

In addition to the changes described above, instrumentation changes can also be performed in this box, although it is recommended that these changes are made with the **Signal Rate** dialog box wherever possible. There are three instrumentation tabs, which are:

Mode: This tab is shown in Figure 6.2 and it permits the user to select the measurement mode, based on the system geometry, for the measurement.

Excitation: This tab is shown in Figure 6.3 and it permits the user to modify the instrumentation associated with the excitation monochromator. The changes that can be made are:

a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.

- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The Light Source radio group is used to select the source to be used in the experiment.
- d. If a computer-controlled iris is fitted, the opening of the iris (as a percentage) can be set using the **Iris Setting** dialog box.
- e. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

Emission: This tab is shown in Figure 6.4 and it permits the user to modify the instrumentation associated with the emission monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The **Detector** radio group is used to select the detector to be used in the experiment.
- d. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

6.2 Card Settings

Depending on the lamp selection, one of the dialog boxes shown in Figure 6.5 and Figure 6.6 will be displayed below the **Manual Measurement** dialog box. If a microsecond flashlamp is used then the measurement will be made with the PCS card operating in MCS (Multi-Channel Scaling) mode and the dialog box shown in Figure 6.5 will be available. The options are:

Time Range: This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 100 μ s to 200 s in sequences of 1-2-4-8.

Channels: This value is the number of channels to be measured over the time range. The available options are 500, 1000, 2000, and 4000 channels.

Lamp Trigger Delay: This value is the delay (in ms) of the lamp trigger with respect to trigger signal to the MCS card that commences an acquisition. It is used to adjust the position the measured data on the data plot so that the data is clearly shown. The setting can be adjusted from 0.001 ms to 90.000 ms in steps of 0.001 ms.

Stop condition: The setting is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Sweeps: This setting stops the measurement after the entered number of sweeps (single measurements) have been completed. Pressing the **Apply** button will display the time required for the entered number of sweeps to be performed.
- **b.** Time: This setting stops the measurement after the specified time has elapsed. Note that the entered time is adjusted by the software to ensure that an integral number of sweeps have been performed. Pressing the **Apply** button will perform this correction without starting the measurement.

c. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.

	Lifetime Sample 1 🛛 🛛 🗙
	Time Range Stop Condition Rates
MCS Measurement 🛛 🔋 🗙	Time Range: 2µs
Time Range Settings	Channels: 1024 💌 Time/ch.: 1.9531 ns
Time Hange: 1s 1.0 ms/ch	
Channels: 1000 💌	Uptions
Lamp Trigger Delay: 5.000 🚔 ms	Time Range Stop Condition Rates
Stop Condition	Time (s) = 10.000 🛓 seconds
Time (s) = 26 🛓 Seconds	🗖 at chan 0 🚊
= 26 Sweeps	
	Time Range Stop Condition Rates
ii	Start Rate: 250 Hz.
	Stop Rate: 260 Hz.
Lamp Trigger Delay: 5.000 ms Stop Condition Time (s) = 26 Seconds = 26 Sweeps Apply	Time Range Stop Condition Rates Time (s) = 10.000 second at chan 0 _ Time Range Stop Condition Rates Start Rate: 250 Hz. Stop Rate: 260 Hz.

Figure 6.5: Multi-Channel Scaling (MCS) measurement options.

Figure 6.6: Time-Correlated Single Photon Counting (TCSPC) options.

If the nanosecond flashlamp or the steady-state lamp are used then measurements will be performed with the TCSPC card and the dialog box shown in Figure 6.6 will be visible. The dialog box has three tabs whose contents are shown in each of the snapshots of the figure. The settings for the **Time Range** and **Stop Condition** are:

Time Range: This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 2.5 ns to 50 μ s in sequences of (after the 2.5 ns value) 1-2-5.

Channels: This value is the number of channels to be measured over the time range. The available options are 512, 1024, 2048, 4096 channels.

Stop condition: The setting is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Time: This setting stops the measurement after the specified time has elapsed.
- b. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.
- c. **Manual:** With this option the user has to stop the measurement manually by pressing the **Stop** button. This will override any entered values of Peak Counts, Time or External.
- d. **External:** This setting requires that an external interrupt signal is required to stop the measurement. This signal is applied to the control port of the TCC-900 card.

The values displayed in the **Rates** tab are read-only with the **Start Rate** value displaying the signal frequency at the Start input of the TCC-900 card and the **Stop Rate** displaying the signal frequency at the Stop input of the TCC-900 card.

Pressing the **Options** button will open a dialog box to change the settings for the TCC-900 card. There are four tabs on the dialog box and these are shown in Figure 6.7.

TCC-900 Options	CFD Time Range Settings Rate Display Mode
CFD Time Range Settings Rate Display Mode Start Stop Threshold: -40 mV Threshold: -300 mV Threshold: -40 mV Zero C.: -10.00 mV Zero C.: -10.00 mV Divider: 1 ▼ Divider: 1	Name: 2µs
OK Cancel Apply	Save Save As Delete
(a) CFD Time Range Settings Rate Display Mode Start Rate Units • Hz • Hz • KHz • MHz • MHz	CFD Time Range Settings Rate Display Mode Acquisition Mode Forward Reverse



CFD

The Constant Fraction discriminator (CFD) set-up is important in detector / electronics optimisation work, in particular when the TCSPC module has been recently installed or when a new detector is to be set up. Both Start and Stop inputs of the TCC-900 card have a CFD to temporally optimise and shape the incoming pulses. Both CFDs can be set up independently. The majority of the CFD parameters can be optimised by software. Only one CFD parameter, the CFD Delay, must be set manually by replacing a delay plug on the TCC900 card. The software settable values are:

a. **Threshold:** The threshold can be set in the range 0 to -2000 mV in steps of 1 mV. As a general rule, select a value about 10% of the average pulse height of the input pulses (pulses from a trigger diode will not have a large pulse height fluctuation, whereas those from a photomultiplier will have a much greater variation).

- b. **Zero Crossing:** The zero crossing can be set in the range of -100 mV to 100 mV in steps of 0.05 mV. As a general rule, start with a small negative value e.g. -3 mV,
- c. Divider: The divider can have values of 1, 2, 4, or 8. Generally, start with a value of 1.

All the above parameters need to be optimised in order to achieve good measurement results, e.g. in respect to the pulse width of the instrument response, to the ratio of the number of dark counts to the number of signal counts, to the elimination of RF and interference problems etc.

The CFD delay also needs to be optimised for the best measurement result. As a general rule the CFD delay should be about 70% of the pulse rise time. Different delay plugs are supplied with the TCC-900 card. The standard CFD delay plugs fitted to both the Start and the Stop channels are 1.0ns and 0.2ns, respectively. If these values are not appropriate for the particular detector, than they can easily be changed. To do this the card needs to be removed from the computer. Once this has been done, the two brown delay plugs (one for the Start the other for the Stop channel) can be located at the edge of the TCC-900 card. To change simply unplug the original plug and insert the new plug.

The **Divider** is not strictly a CFD parameter, but it may be needed when the TCC-900 card is operating in **Reverse** mode. By using the original (laser) repetition rate in reverse TCSPC operation only one period of the high repetition signal can be seen. This may be inconvenient when measuring in a time range close to the inverse of the (laser) repetition rate. By changing the divider to either 2, 4, or 8, more than just one signal periods can be measured. Note that the settings of the Divider will not have an impact on the displayed signal rate, it will only be apparent when acquiring the lifetime data.

Time Range Settings

When the TCC-900 card has been set for the first time and one is at the stage that signal rates are displayed, the most likely difficulty is firstly finding the measurement on the measurement display and secondly positioning the rising edge close to the left side of the window. Changing either of the two required shift parameters will shift the entire measurement (decay and/or IRF measurement) either to the left or to the right on the measurement window. The parameters are:

- a. Shifting Delay: This is essentially a series of wire delays integrated in the TCC-900 card. Generally, for large time ranges (above 100 ns) a shifting delay is required in order to shift the measurement to the right to cover the full lifetime measurement, including the rising edge. For short time ranges a small shifting delay may be required in order to compensate for the different cable lengths in the start and stop channels. By correctly "placing" the measurement, full use of the x-axis resolution can be made. A large shifting delay is further required when working in Reverse mode. In this mode the exciting light pulse (usually a laser pulse) needs to be shifted in time to arrive at the TC-900 card later than latest expected decay photon, i.e. the shifting delay must be slightly longer than the selected time range.
- b. **TAC Offset:** The TAC offset can also be used to shift the measurements. The default setting is 500 ns. Values between 300 ns and 1000 ns should only be used for time ranges above 5ns. For the two shortest time ranges (2.5 ns and 5.0 ns) the offset spans from 1200 ns to 4000 ns.

Once the **Time Range Settings** box has been closed the selected delay value will always be active for the particular time range. Other time ranges maintain the old delay values.

Rate Display

The **Rate Display** tab permits the changing of the format of the count rate (per seconds) on the main data acquisition screen. The frequency counters on the TCC-900 card can not measure frequencies above 10 MHz correctly. When working with high synchronisation frequencies either kHz or MHz should be chosen, otherwise the display will be wrong.

Mode

This tab allows the user to change between **Forward** and **Reverse** mode. The change will affect the x-axes of the display, which is effectively reversed when changing between the two modes. If one chooses to work in the Reverse mode, the laser synchronisation frequency (the higher of the two rates) must be connected to the Stop input of the TCC-900 card. Alternatively when working in the forward mode the source synchronisation frequency needs to be connected to the Start channel.

6.3 Multiple

It is possible to perform a sequence of measurements within a single sequence and the multiple lifetime measurement dialog box (shown in Figure 6.8) is used for this purpose.

Multiple Lifetime Measurem 🙎	×
Set up Sample Position Measure Samples Image: Contract of the set of the	
✓ Interleave sample	es
Sample Set Up Time Set Up	
Instrument Response	
Measure this sample 5 📑 times	
🔽 Save data each time	
Wait for 0.0 🚔 minutes after each cycle	
Start Cance	1

Figure 6.8: Dialog box for multiple lifetime measurements.

The constituents of this dialog box are:

Set up Sample Position: This radio group is use to select the sample position that is to be modified by pressing the Sample Set Up button. The options are only available if the multiple sample holder has been installed.

Measure Samples: To include the sample in a particular position, the corresponding check box should be clicked.

Interleave samples: Interleaving affects the order in which samples are measured during automated lifetime measurements. If the measurements are interleaved, the first measurement is made from each of the selected samples, then the second from each, then the third from each, and so on until all acquisitions are made from all samples. If one of the samples is specified as an instrument response, acquisitions from this container will be copied into the containers of the sample measurements on completion. If the measurements are not interleaved, all measurements are completed with a selected sample before passing on to the next. There is no combination of instrument response data with other containers.

Instrument Response: Checking this box will indicate to the software that the measurement sequence is for an instrument response scan and the plotted data will be flagged accordingly.

Measure this sample...: This edit box permits the user to enter the number of times that a particular sample should be measured.

Save data each time: Checking this box will save data each time a measurement is made and will display the measurements separately. Alternatively, the measurements will be aggregated together into a single plot.

Wait for...: Entering a value in this box will pause the measurement for the specified time. This is to permit the user to make changes to the sample(s), for example in chemical concentrations.

Pressing the **Start** button will commence the measurement and pressing the **Cancel** button will close the dialog box without performing any measurements.

6.3.1 Sample Set Up Button

Pressing this button will open the dialog box shown in Figure 6.9.

Sample Set up 🛛 📪 🔀	Mode Excitation Emission 1
Set up Sample © 1 C 2 C 3	Wavelength (nm) Increment (nm) 300.00 — 0.05 ▼
Mode Excitation Emission 1 Measurement Mode • L1 C L2 C T	Angle: 0.0 V
	Iris Setting: 5 🚖 Δλ: 0.18 nm
	(b)
	Mode Excitation Emission 1
	Mode Excitation Emission 1 Wavelength (nm) Increment (nm)
OK Cancel	Mode Excitation Emission 1 Wavelength (nm) Increment (nm) 400.00 - 0.05 -
OK Cancel	Mode Excitation Emission 1 Wavelength (nm) Increment (nm) 400.00 Polariser Detector Polariser Detector
(a)	Mode Excitation Emission 1 Wavelength (nm) Increment (nm) 400.00 + 0.05 + Polariser Detector Angle: 0.0 + PCS 1
(a)	Mode Excitation Emission 1 Wavelength (nm) Increment (nm) 400.00 ♀ 0.05 ▼ Polariser Detector Angle: 0.0 ▼ △X: 1.50

Figure 6.9: Multiple measurement sample set up dialogs with (a) Mode tab, (b) Excitation tab and (c) Emission tab shown.

There are three instrumentation tabs, which are:

Mode: This tab is shown in Figure 6.9(a) and it permits the user to select the measurement mode, based on the system geometry, for the measurement.

Excitation: This tab is shown in Figure 6.9(b) and it permits the user to modify the instrumentation associated with the excitation monochromator. The changes that can be made are:

- f. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- g. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- h. The Light Source radio group is used to select the source to be used in the experiment.
- i. If a computer-controlled iris is fitted, the opening of the iris (as a percentage) can be set using the **Iris Setting** dialog box.
- j. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

Emission: This tab is shown in Figure 6.9(c) and it permits the user to modify the instrumentation associated with the emission monochromator. The changes that can be made are:

- e. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- f. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- g. The Detector radio group is used to select the detector to be used in the experiment.
- h. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

6.3.2 Time Set Up Button

Pressing the **Time Set Up button** opens either the dialog box shown in Figure 6.10 (for measurements made with the steady-state or nanosecond flashlamps) or Figure 6.11 (for measurements made with the microsecond flashlamp).

The settings for the Time Range and Stop Condition tabs are:

Time Range: This tab is used to set the timing conditions for the measurement.

- a. **Time Range:** This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 2.5 ns to 50 μ s in sequences of (after the 2.5 ns value) 1-2-5.
- b. **Channels:** This value is the number of channels to be measured over the time range. The available options are 512, 1024, 2048, 4096 channels.

Stop condition: The tab is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Time: This setting stops the measurement after the specified time has elapsed.
- b. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.

- c. **Manual:** With this option the user has to stop the measurement manually by pressing the **Stop** button. This will override any entered values of Peak Counts, Time or External.
- d. **External:** This setting requires that an external interrupt signal is required to stop the measurement. This signal is applied to the control port of the TCC-900 card.

Lifetime Sample 1 🛛 🗙	Time Range Stop Condition
Time Range Stop Condition	Peak Counts 👻 = 10000 🏯 counts
Time Range: 2µs	T at chan
Channels: 1024 💌 Time/ch.: 1.9531 ns	
Options OK Cancel	
(a)	(b)

Figure 6.10: Multiple measurement time set up dialogs for TCSPC mode measurements showing (a) Time Range tab and (b) Stop Condition tab.

Pressing the Options button will open the TCC-900 settings dialog box. This is described in Section 6.2.

MCS Measurement 💦 👔
Time Range Settings
Time Range: 400ms 💌
Channels: 1000 💌
Lamp Trigger Delay: 0.100 📩 ms
Stop Condition
Time (s) 💌 = 10 🐳 Seconds
= 10 Sweeps
OK Cancel

Figure 6.11: Multiple measurement time set up dialog box for MCS measurements.

The items in the MCS Measurement dialog box are:

Time Range: This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 100 μ s to 200 s in sequences of 1-2-4-8.

Channels: This value is the number of channels to be measured over the time range. The available options are 500, 1000, 2000, and 4000 channels.

Lamp Trigger Delay: This value is the delay (in ms) of the lamp trigger with respect to trigger signal to the MCS card that commences an acquisition. It is used to adjust the position the measured data on the data plot

so that the data is clearly shown. The setting can be adjusted from 0.001 ms to 90.000 ms in steps of 0.001 ms.

Stop condition: The setting is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Sweeps: This setting stops the measurement after the entered number of sweeps (single measurements) have been completed. Pressing the Apply button will display the time required for the entered number of sweeps to be performed.
- **b.** Time: This setting stops the measurement after the specified time has elapsed. Note that the entered time is adjusted by the software to ensure that an integral number of sweeps have been performed. Pressing the **Apply** button will perform this correction without starting the measurement.
- c. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.

6.4 TRES

Time-Resolved Emission Scans (TRES) are used to examine how lifetime decay change with wavelength. They are performed by measuring decay over a pre-defined set of equally-spaced wavelengths. The dialog boxes use for this purpose are shown in Figure 6.12.

TRES Excitation Measureme ? × Sample Measure Samples 1 2 3 Image: The sample sampl	Sample Measurements No Image: 1 1 2 3 Image: 1 1 2 3 Image: 1 1 1 1 Image: 1 1 1 1 1 Image: 1 1 1 1 1 1 Image: 1 1 1 1 1 1 1 Image: 1 1 1 1 1 1 1 1 1 Image: 1 1
Sample Set Up Time Set Up Instrument Response Wavelength Setup (nm) Ex 10000 + to 420.00 + Step 10.00 + Em1 200.00 +	Sample Set Up Instrument Response Wavelength Setup (nm) Ex SSSS Em1 592.00 to 593.00 Step 0.05
Start Cancel	Start Cancel
(a)	(b)

Figure 6.12: Dialog boxes for measuring (a) TRES excitation scans and (b) TRES emission scans.

The common constituents of these dialog boxes are:

Set up Sample Position: This radio group is use to select the sample position that is to be modified by pressing the **Sample Set Up** button. The options are only available if the multiple sample holder has been installed.

Measure Samples: To include the sample in a particular position, the corresponding check box should be clicked.

Interleave samples: Interleaving affects the order in which samples are measured during automated lifetime measurements. If the measurements are interleaved, the first measurement is made from each of the selected samples, then the second from each, then the third from each, and so on until all acquisitions are made from all samples. If one of the samples is specified as an instrument response, acquisitions from this container will be copied into the containers of the sample measurements on completion. If the measurements are not interleaved, all measurements are completed with a selected sample before passing on to the next. There is no combination of instrument response data with other containers.

Instrument Response: Checking this box will indicate to the software that the measurement sequence is for an instrument response scan and the plotted data will be flagged accordingly.

Wavelength Setup: For an Excitation TRES measurement the wavelength of the emission monochromator is fixed whilst the wavelength of the excitation monochromator is varied. The start, stop and step (increment) wavelengths are entered in this group box, as can be seen in Figure 6.12(a). For emission TRES measurements it is the excitation monochromator wavelength that is allowed to change and the emission monochromator whose wavelengths are changed. The box to do this is shown in Figure 6.12(b).

Pressing the **Start** button will commence the measurement and pressing the **Cancel** button will close the dialog box without performing any measurements.

6.4.1 Sample Set Up Button

Pressing this button will open the dialog box shown in Figure 6.13.

There are three instrumentation tabs, which are:

Mode: This tab is shown in Figure 6.13(a) and it permits the user to select the measurement mode, based on the system geometry, for the measurement.

Excitation: This tab is shown in Figure 6.13(b) and it permits the user to modify the instrumentation associated with the excitation monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The Light Source radio group is used to select the source to be used in the experiment.
- d. If a computer-controlled iris is fitted, the opening of the iris (as a percentage) can be set using the **Iris Setting** dialog box.
- e. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

Sample Set up 🛛 📪 🗙	Mode Excitation Emission 1
Set up Sample © 1 © 2 © 3	Wavelength (nm) Increment (nm) 300.00 🐳 0.05 💌
Mode Excitation Emission 1 Measurement Mode • L1 • L2 • T	Polariser Light Source Angle: 0.0 Tris Setting: 5 Δλ: 0.18
	(b) Mode Excitation Emission 1
OK Cancel	Wavelength (nm) Increment (nm) 400.00
(a)	Angle: 0.0 C PCS 1 Δλ: 1.50 nm
	(c)

Figure 6.13: Multiple measurement sample set up dialogs with (a) Mode tab, (b) Excitation tab and (c) Emission tab shown.

Emission: This tab is shown in Figure 6.13(c) and it permits the user to modify the instrumentation associated with the emission monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The Detector radio group is used to select the detector to be used in the experiment.
- d. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

6.4.2 Time Set Up Button

Pressing the **Time Set Up button** opens either the dialog box shown in Figure 6.14 (for measurements made with the steady-state or nanosecond flashlamps) or Figure 6.23 (for measurements made with the microsecond flashlamp).

Lifetime Sample 1 🛛 🛛 🗙	Time Range Stop Condition
Time Range Stop Condition	Peak Counts 💌 = 10000 🛖 counts
Time Range: 2µs	at chan 0
Channels: 1024 💌 Time/ch.: 1.9531 ns	
Options OK Cancel	
(a)	(b)

Figure 6.14: Multiple measurement time set up dialogs for TCSPC mode measurements showing (a) Time Range tab and (b) Stop Condition tab.

The settings for the Time Range and Stop Condition tabs are:

Time Range: This tab is used to set the timing conditions for the measurement.

- a. **Time Range:** This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 2.5 ns to 50 μ s in sequences of (after the 2.5 ns value) 1-2-5.
- b. **Channels:** This value is the number of channels to be measured over the time range. The available options are 512, 1024, 2048, 4096 channels.

Stop condition: The tab is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Time: This setting stops the measurement after the specified time has elapsed.
- b. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.
- c. **Manual:** With this option the user has to stop the measurement manually by pressing the **Stop** button. This will override any entered values of Peak Counts, Time or External.
- d. **External:** This setting requires that an external interrupt signal is required to stop the measurement. This signal is applied to the control port of the TCC-900 card.

Pressing the **Options** button will open the TCC-900 settings dialog box. This is described in Section 6.2.

The items in the MCS Measurement dialog box are:

Time Range: This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 100 μ s to 200 s in sequences of 1-2-4-8.

Channels: This value is the number of channels to be measured over the time range. The available options are 500, 1000, 2000, and 4000 channels.

Lamp Trigger Delay: This value is the delay (in ms) of the lamp trigger with respect to trigger signal to the MCS card that commences an acquisition. It is used to adjust the position the measured data on the data plot so that the data is clearly shown. The setting can be adjusted from 0.001 ms to 90.000 ms in steps of 0.001 ms.

MCS Measurement	X
Time Range Settings	
Time Range: 400ms 💌	leh
Channels: 1000 💌	Ch
Lamp Trigger Delay: 0.100 📑 ms	
Stop Condition	
Time (s) 🗨 = 10 🜩 Secor	nds
= 10 Sweeps	
OK Cance	

Figure 6.15: Multiple measurement time set up dialog box for MCS measurements.

Stop condition: The setting is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- **a. Sweeps:** This setting stops the measurement after the entered number of sweeps (single measurements) have been completed. Pressing the **Apply** button will display the time required for the entered number of sweeps to be performed.
- **b.** Time: This setting stops the measurement after the specified time has elapsed. Note that the entered time is adjusted by the software to ensure that an integral number of sweeps have been performed. Pressing the **Apply** button will perform this correction without starting the measurement.
- c. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.

6.5 Anisotropy Lifetime

Anisotropy lifetime measurements are used to examine the temporal development of fluorescence anisotropy. The dialog box used for this purpose is shown in Figure 6.16.

The common constituents of these dialog boxes are:

Sample: This radio group is use to select the sample position that is to be modified by pressing the **Sample Set Up** button. The options are only available if the multiple sample holder has been installed.

Instrument Response: Checking this box will indicate to the software that the measurement sequence is for an instrument response scan and the plotted data will be flagged accordingly.

Measure this Sample: Enter the number of times that the sample has to be measured.

Wait for: Enter the interval between consecutive measurements. This option allows the user to make changes to the sample between measurements.

Polariser Repetition: Enter the number of polariser repetition sequences in this edit box.

Anisotropy Lifetime Measurements ? × Sample © 1 © 2 © 3
Sample Set Up Time Set Up
Instrument Response
Measure this Sample 1 🚆 times
Wait for 0.00 🔺 minutes between cycles
Polariser Repetition 1 🚔
Polariser Sequence
C ExV, Em1 V, Em2 H
C ExV, Emily then H, Em21(then H
G Factor
Measure now by integrating TCSPC
○ Measure from signal rate for 1 seconds
○ Fixed value of 0.00 🚔
Cancel Start

Figure 6.16: Dialog box for measuring anisotropy lifetime scans.

Polariser sequence: If the vertical polarisation direction is selected for the excitation channel the spectrally resolved fluorescence intensities I_{VV} and I_{VH} will be sequentially measured. If the horizontal polarisation direction has been selected the excitation scans I_{HV} and I_{HH} , required for G-factor correction, will be sequentially measured. In **T-geometry**, six different modes of automatic polarised fluorescence measurements can be set up, with three of them used to measure curves in order to determine the G-factor.

If the vertical polarisation direction is selected for the excitation channel and the first of the three options is selected for the emission channels then I_{VV} and I_{VH} will be respectively and simultaneously measured at the two emission arms.

If the vertical polarisation direction is selected for the excitation channel and the second of the 3 options is selected for the emission channels I_{VV} and I_{VH} will be measured at emission arm 1 subsequently, whereas a user defined polarised measurement will be made simultaneously in the second emission arm. If the vertical polarisation direction is selected for the excitation channel and the last of the 3 options is selected for the emission channels the spectrally resolved fluorescence intensities I_{VV} will be measured at both emission arms simultaneously and I_{VH} will be measured for both arms afterwards.

G Factor: The measurements for horizontal oriented polariser in the excitation channel are similar to the measurements described above. They are used to make the G factor correction..

All polarised measurements are stored automatically with the file extensions ".VV", ".VH", ".HV", and ".HH", dependent upon the orientation of the excitation and emission polarisers.
Pressing the **Start** button will commence the measurement and pressing the **Cancel** button will close the dialog box without performing any measurements.

6.5.1 Sample Set Up Button

Pressing this button will open the dialog box shown in Figure 6.17.

There are three instrumentation tabs, which are:

Mode: This tab is shown in Figure 6.9(a) and it permits the user to select the measurement mode, based on the system geometry, for the measurement.

Excitation: This tab is shown in Figure 6.17(b) and it permits the user to modify the instrumentation associated with the excitation monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The Light Source radio group is used to select the source to be used in the experiment.
- d. If a computer-controlled iris is fitted, the opening of the iris (as a percentage) can be set using the **Iris Setting** dialog box.
- e. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

Emission: This tab is shown in Figure 6.17(c) and it permits the user to modify the instrumentation associated with the emission monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The **Detector** radio group is used to select the detector to be used in the experiment.
- d. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

Sample Set up 🛛 🔋 🗙	Mode Excitation Emission 1
Set up Sample © 1 C 2 C 3	Wavelength (nm) Increment (nm) 300.00 ⊕ 0.05 ▼
Mode Excitation Emission 1 Measurement Mode © L1 C L2 C T	Angle: 0.0 V
	Iris Setting: <mark>5 🚔 Δλ: 0.18 nm</mark>
	(b)
	Mode Excitation Emission 1
	Wavelength (nm) Increment (nm)
OK Cancel	400.00 🛫 0.05 💌
	Polariser Detector
(a)	Angle: 0.0 C PCS 1
	Δλ: 1.50 nm
	(c)

Figure 6.17: Multiple measurement sample set up dialogs with (a) Mode tab, (b) Excitation tab and (c) Emission tab shown.

6.5.2 Time Set Up Button

Pressing the **Time Set Up button** opens either the dialog box shown in Figure 6.18 (for measurements made with the steady-state or nanosecond flashlamps) or Figure 6.19 (for measurements made with the microsecond flashlamp).

Lifetime Sample 1 🛛 🗙	Time Range Stop Condition
Time Range Stop Condition	Peak Counts 💌 = 10000 📥 counts
Time Range: 2µs	T at chan 0
Channels: 1024 💌 Time/ch.: 1.9531 ns	
Options OK Cancel	
(a)	(b)

Figure 6.18: Multiple measurement time set up dialogs for TCSPC mode measurements showing (a) Time Range tab and (b) Stop Condition tab.

The settings for the Time Range and Stop Condition tabs are:

Time Range: This tab is used to set the timing conditions for the measurement.

- a. **Time Range:** This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 2.5 ns to 50 μ s in sequences of (after the 2.5 ns value) 1-2-5.
- b. **Channels:** This value is the number of channels to be measured over the time range. The available options are 512, 1024, 2048, 4096 channels.

Stop condition: The tab is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Time: This setting stops the measurement after the specified time has elapsed.
- b. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.
- c. **Manual:** With this option the user has to stop the measurement manually by pressing the **Stop** button. This will override any entered values of Peak Counts, Time or External.
- d. **External:** This setting requires that an external interrupt signal is required to stop the measurement. This signal is applied to the control port of the TCC-900 card.

Pressing the Options button will open the TCC-900 settings dialog box. This is described in Section 6.2.

MCS Measurement 🛛 🙎 🗙
Time Range Settings
Time Range: 400ms 💌 0.4 ms/ch
Channels: 1000 💌
Lamp Trigger Delay: 0.100 🛖 ms
Stop Condition
Time (s) 💌 = 10 🛨 Seconds
= 10 Sweeps
OK Cancel

Figure 6.19: Multiple measurement time set up dialog box for MCS measurements.

The items in the MCS Measurement dialog box are:

Time Range: This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 100 μ s to 200 s in sequences of 1-2-4-8.

Channels: This value is the number of channels to be measured over the time range. The available options are 500, 1000, 2000, and 4000 channels.

Lamp Trigger Delay: This value is the delay (in ms) of the lamp trigger with respect to trigger signal to the MCS card that commences an acquisition. It is used to adjust the position the measured data on the data plot so that the data is clearly shown. The setting can be adjusted from 0.001 ms to 90.000 ms in steps of 0.001 ms.

Stop condition: The setting is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Sweeps: This setting stops the measurement after the entered number of sweeps (single measurements) have been completed. Pressing the **Apply** button will display the time required for the entered number of sweeps to be performed.
- **b.** Time: This setting stops the measurement after the specified time has elapsed. Note that the entered time is adjusted by the software to ensure that an integral number of sweeps have been performed. Pressing the **Apply** button will perform this correction without starting the measurement.
- c. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.

6.6 Temperature Lifetime

Temperature lifetime scans are used to examine the effect of temperature on temporal development of a fluorescence decay. The dialog boxes use for this purpose are shown in Figure 6.20.

Temperature Lifetime Meas ? 🗙
Sample © 1 © 2 © 3
Sample Set Up
Instrument Response
Temperature (K) Setup
From 295.00 🛨 to 300.00 🛨 step 1.00 🛨
Accuracy Band (±) 0.10 🚔 🔽 Auto PID
Stabilisation Time 0 🚖 s
Start Cancel

Figure 6.20: Dialog box for measuring temperature lifetime scans.

The common constituents of these dialog boxes are:

Sample: This radio group is use to select the sample position that is to be modified by pressing the **Sample Set Up** button. The options are only available if the multiple sample holder has been installed.

Instrument Response: Checking this box will indicate to the software that the measurement sequence is for an instrument response scan and the plotted data will be flagged accordingly.

Scan From/To/Step: Enter the temperature range of interest in Kelvin (Scan from/to) and the number of degrees Kelvin between each measurement (Step).

Accuracy Band: Enter the temperature tolerance range (in Kelvin) within which is the temperature is assumed to be stable.

Stabilisation Time: Enter the time that the cryostat has to stay at each temperature before a measurement is acquired.

Auto PID: The Proportional, Integral and Derivative settings can either be set manually on the cryostat controller or selected automatically from the learnt characteristics of the system by means of the **Auto PID** check box in the **Temperature Scan Parameters**. When the cryostat temperature is set to go to a higher temperature, the temperature increase has a hysteresis profile. This is one of the reasons for a stabilisation time. In order to minimise the hysteresis while maximising the speed of the temperature increase, the proportional, integral and derivative settings can be altered. We recommend that the temperature controller automatically sets these. More information on this can be found in the cryostat instruction manuals.

Pressing the **Start** button will commence the measurement and pressing the **Cancel** button will close the dialog box without performing any measurements.

6.6.1 Sample Set Up Button

Pressing this button will open the dialog box shown in Figure 6.21.

There are three instrumentation tabs, which are:

Mode: This tab is shown in Figure 6.21(a) and it permits the user to select the measurement mode, based on the system geometry, for the measurement.

Excitation: This tab is shown in Figure 6.21(b) and it permits the user to modify the instrumentation associated with the excitation monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The Light Source radio group is used to select the source to be used in the experiment.
- d. If a computer-controlled iris is fitted, the opening of the iris (as a percentage) can be set using the **Iris Setting** dialog box.
- e. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

Sample Set up 🛛 🔋 🗙	Mode Excitation Emission 1
Set up Sample © 1 O 2 O 3	Wavelength (nm) Increment (nm) 300.00 🐳 0.05 💌
Mode Excitation Emission 1 Measurement Mode © L1 © L2 © T	Angle: 0.0 V 0.10
	(b)
OK Cancel	Wavelength (nm) Increment (nm)
(a)	Angle: 0.0 C PCS 1
	∆: 1.50 nm (c)

Figure 6.21: Multiple measurement sample set up dialogs with (a) Mode tab, (b) Excitation tab and (c) Emission tab shown.

Emission: This tab is shown in Figure 6.21(c) and it permits the user to modify the instrumentation associated with the emission monochromator. The changes that can be made are:

- a. The monochromator wavelength can be changed in the **Wavelength** edit box by either entering a new wavelength directly into the box or by clicking on the increment arrows. The size of the increment is determined by the value of the **Increment** drop-down box.
- b. If fitted, the polariser setting can be changed by clicking the **Angle** drop-down box and choosing the desired angle from the options available.
- c. The Detector radio group is used to select the detector to be used in the experiment.
- d. If computer-controlled slits have been fitted then their separation is sown in the bottom, righthand corner of the tab, although the value cannot be changed here. This can only be done in the **Signal Rate** dialog box.

6.6.2 Time Set Up Button

Pressing the **Time Set Up button** opens either the dialog box shown in Figure 6.22 (for measurements made with the steady-state or nanosecond flashlamps) or Figure 6.23 (for measurements made with the microsecond flashlamp).

Lifetime Sample 1 🛛 🛛 🗙	Time Range Stop Condition
Time Range Stop Condition	Peak Counts 💌 = 10000 🛖 counts
Time Range: 2µs	at chan 0
Channels: 1024 💌 Time/ch.: 1.9531 ns	
Options OK Cancel	
(a)	(b)

Figure 6.22: Multiple measurement time set up dialogs for TCSPC mode measurements showing (a) Time Range tab and (b) Stop Condition tab.

The settings for the Time Range and Stop Condition tabs are:

Time Range: This tab is used to set the timing conditions for the measurement.

- a. **Time Range:** This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 2.5 ns to 50 μ s in sequences of (after the 2.5 ns value) 1-2-5.
- b. **Channels:** This value is the number of channels to be measured over the time range. The available options are 512, 1024, 2048, 4096 channels.

Stop condition: The tab is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Time: This setting stops the measurement after the specified time has elapsed.
- b. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.
- c. **Manual:** With this option the user has to stop the measurement manually by pressing the **Stop** button. This will override any entered values of Peak Counts, Time or External.
- d. **External:** This setting requires that an external interrupt signal is required to stop the measurement. This signal is applied to the control port of the TCC-900 card.

Pressing the Options button will open the TCC-900 settings dialog box. This is described in Section 6.2.

MCS Measurement	? ×
Time Range Settings	
Time Range: 400ms 💌	4 ms/ch
Channels: 1000 💌	4 mar on
Lamp Trigger Delay: 0.100 🚔 ms	
Stop Condition	
Time (s) 💌 = 10 🚔 9	Seconds
= 10 Sweeps	
ОК	Cancel

Figure 6.23: Multiple measurement time set up dialog box for MCS measurements.

The items in the MCS Measurement dialog box are:

Time Range: This is the overall duration of a single measurement and is selectable from the options in the drop-down box. The values range from 100 μ s to 200 s in sequences of 1-2-4-8.

Channels: This value is the number of channels to be measured over the time range. The available options are 500, 1000, 2000, and 4000 channels.

Lamp Trigger Delay: This value is the delay (in ms) of the lamp trigger with respect to trigger signal to the MCS card that commences an acquisition. It is used to adjust the position the measured data on the data plot so that the data is clearly shown. The setting can be adjusted from 0.001 ms to 90.000 ms in steps of 0.001 ms.

Stop condition: The setting is used to determine the stop condition for the measurements. There are three possible ways of setting the stop condition and these are:

- a. Sweeps: This setting stops the measurement after the entered number of sweeps (single measurements) have been completed. Pressing the **Apply** button will display the time required for the entered number of sweeps to be performed.
- b. Time: This setting stops the measurement after the specified time has elapsed. Note that the entered time is adjusted by the software to ensure that an integral number of sweeps have been performed. Pressing the Apply button will perform this correction without starting the measurement.
- c. **Peak Counts:** The setting permits the user to stop the measurement once a specified number of counts have been received on the (by default) highest amplitude channel of the data. If desired, the channel that the peak counts are measured can be specified by the user.

7 Viewing Data

The **View** menu (described in the next chapter) offers options for displaying the data. This menu will only be available if data has previously been loaded from disk or if new data has been acquired. It is possible to view the acquired data in up to four different ways. These are:

- a. As 2D plot
- b. As 3D plot
- c. As Text plot
- d. As Contour plot

The **As 2D Plot** and **As Text** options will always be available but **As 3D Plot** and **As Contour Plot** are only available if the active window contains a map (emission or absorption) measurement.

The Plot Options depend on the type of plot currently active and are used to change the appearance of a plot. To display the plot options menu go to the View Menu and click in the "Plot Options menu". You can also display this menu by clicking the **Plot Options** icon in the toolbar.

7.1 As 2D plot

This is the default method and is also the only one that you can be used during data acquisition. It provides a two dimensional plot of the data being acquired. It also provides unique features such as a Cursor and Peak Searching. This 2D plot representation is also available by clicking the **View as 2D Plot** icon button. An example of a 2D plot is shown in Figure 7.1.



Figure 7.1: Example of a 2-D plot (emission scan of Anthracene).

Pressing the **Plot Options** menu item or toolbar button when a 2D plot is being displayed will open the dialog shown in Figure 7.2.

Current 2D Plot Option	ns 🛛 🗙
General Lifetime	
Font Arial	Change
Screen Colour Options	
Background 💌	-Miscellaneous Options
Colour	ge Show Grid
Line Style Line, Solid	Y
Line Width 1	
Save as <u>D</u> efault OK	<u>C</u> ancel Apply

Figure 7.2: General 2D Plot menu

The changes that can be made in this menu are:

Font: Click the **Change** button inside the Font box. When the dialog box opens, select the font style and size as desired.

Screen Colour Options: This submenu allows you to select the part of the plot you want to change in the item list (background, traces, axis ,etc). From the drop-down box select the item whose coulour is to be changed and then press the **Change** button to select a new colour.

Line Style: Select a new line style from the options available.

Line Width: Enter a new line width in the edit box, either directly, or by using the arrows at the side of the box.

Miscellaneous Options: The **Show Grid** checkbox determines whether the plot should have a grid superimposed on it and the **Show Cursor** checkbox determines whether a cursor should be visible on the plot.

To activate any changes, the **Apply** button should be pressed. After a change has been made, the user can select to make these changes as the default option for plotting. To do this, click the **Save as Default** button. For the case when no measurements are open and it is desired to set the defaults, it can be done in the same way by going into **Plot Defaults** in the **Options** menu.

Depending on the 2D-plot type, the second tab in the **Plot options menu** will display a tab menu according to the plot type. Thus, there can be spectral, lifetime and fit tab menus. These are shown in Figure 7.3.

In these tab menus, the user can select and modify the X and Y-scaling modes, and changes to the axis label.

X-axis: For a spectral plot type, the only X-axis available is Wavelength. For a lifetime or fit type, it is possible to select Time or Channels as the Y-axis.

Y-axis: In all the 2D-plot types, the Y-axis can be changed from linear to logarithmic, although the logarithmic scale may not be possible depending on the Y-range. The same effect as switching between linear and logarithmic Y-scale for an active window can be obtained by clicking the icons called "lin" and "log" in the toolbar.

Scale from Zero: If this option is selected the Y-axis range will begin at zero.



Figure 7.3: Tab menus for (a) Lifetime 2-D plot , (b) Spectral 2-D plot, and (c) Fit 2-D plot.

7.2 As 3D plot

This option allows a measurement containing several scans to be viewed in three dimensions. The data can be viewed from various angles and heights to get a fuller picture of the measurement results. Any measurement can be viewed as a 3D plot as long as it has the properties of a map i.e. it contains more than one scan and all scans have equal step sizes.



Figure 7.4: Example of a 3D representation of a map measurement.

If the active window is a 3D plot, the plot options available can be summarised in the following list:

- a. Viewing Position
- b. Appearance
- c. Data
- d. Axes
- e. Printing

Viewing position: This allows the user to select the angle (**View From**) and height (**View Height**) from which to view the data. The centre position provides an overhead view, where only the **High** viewing height is used. All viewing positions look towards the centre of the data.

3D Plot Options	? ×
Viewing Position Appearance	Data Axes Printing
View from	View Height
Back	 High
Left C C Right	C Middle
Front	C Base
Make Default OK	Cancel Apply

Figure 7.5: Viewing Position tab.



Figure 7.6: Two views of a 3D plot. (a) View from front-right angle and middle height and (b) from front angle and middle height .

Appearance: In this menu the user can switch between perspective and orthogonal viewing modes. The **Keep** Aspect option is always on in **Perspective** mode. Here, you can also select the colours for each of the components of the plot (background, axes, etc) and the font type and size. To change the lighting, select **Light1** or **Light2** in the colours list. On the **Colour** dialog, select the colour you want for that light. To change the luminosity (or brightness) of the light, use the slider on the right hand side of the Colour dialog. **Light 1** is positioned in line with the left axis wall, in front of the x-scale. **Light 2** is positioned in line with the back axis wall, to the right of the z-scale.

3D Plot Options ? 🗙	3D Plot Options
Viewing Position Appearance Data Axes Printing	Viewing Position Appearance Data Axes Printing
Perspective Keep aspect ratio Colours Background Font Arial Change	 Draw Contours Style Filled Surface Triangle Mesh Dots
Make Default OK Cancel Apply	Make Default OK Cancel Apply
(a)	(b)

Figure 7.7: Appearance (a) and Data (b) tabs.

Data: This tab menu permits the choice the surface style for the data. The options are a filled surface, mesh surface or dots.

Axes: Tick the box labelled Filled Axis Panels for solid axis panels, clear it for lines.

Viewing Position Appearance Data Axes Printing Viewing Image: Filled Axis Panels Filled Axis Panels Print Filled Axis Panels Print Filled Axis Panels	Position Appearance Data Axes Printing esolution Default Printer: Panasonic KX-P6300 GDI					
✓ Filled Axis Panels ✓ Filled Axis Panels ✓ Show Map Container's Z Axis as Description Labels	esolution Default Printer: Panasonic KX-P6300 GDI					
Show Map Container's Z Axis as Description Labels	Default Printer: Panasonic KX-P6300 GDI					
-Y-Axis	Default Printer: Panasonic KX-P6300 GDI Page Resolution: 2716 x 3917 pixels					
C Logarithmic						
Si	eed Quality					
Make Default OK Cancel Apply Make Default	fault OK Cancel Apply					
(a)	(h)					

Figure 7.8: Axes (a) and Printing (b) tabs

Printing: On the **Printing** tab, slide the pointer until the desired print resolution is displayed. The 3D plot will always appear the same size on the printed page regardless of this setting. Sliding the pointer fully to the right means that the 3D plot will be printed at the printer maximum resolution. This results in the creation of a very large 3D-plot bitmap, which requires a lot of memory. If this causes Windows to invoke virtual memory on disk, printing may take several minutes. Printing times may be reduced by decreasing the print resolution, using 256-colour mode instead of 24-bit true colour, or by adding more memory to the system. The slider position second from the left is the default.

7.3 As Text

Selecting this option displays the measurement data as a table of values and allows each of them to be edited individually if desired. A plot in text mode is shown on in Figure 7.9; the upper part displays measurement conditions, i.e. scan properties; while the lower part, a table, displays the measurement data. The data can be modified only in this mode. If it is wished to modify a datum value, click the corresponding cell in the table and type a new value. The plot options for Text mode are the same as those for 2D.

Scan : Em Wave	elength:	Em 368	Ant ission Scan 3.00 to 500.0	thracene EN Experi : M300 00 step 1.00	/ M300_1(E i mental Cond) nm	mission Sca ditions	n)				
Ex Wave Em Slit S Em Polar Dwell Tin	Length: Size: riser: ne:	358.00 nm 1.00 mm Ex Slit Size: No polariser Ex Polariser: 0.20 S Temperature:					0.50 mm No polariser				
Data / 10e3 (Counts)											
nm	0	1	2	3	4	5	6	7	8	9	
368	0.133	0.156	0.199	0.279	0.409	0.725	1.210	1.879	2.480	2.815	
378	2.739	2.320	2.031	1.888	1.778	1.610	1.380	1.138	0.941	0.873	
388	0.737	0.763	0.779	0.855	1.032	1.339	1.694	2.093	2.573	3.005	
398	3.221	3.354	3.144	2.882	2.659	2.192	1.917	1.809	1.491	1.354	
408	1.222	1.004	0.916	0.832	0.760	0.780	0.802	0.872	0.982	1.050	
418	1.274	1.388	1.504	1.594	1.661	1.621	1.666	1.503	1.361	1.261	
428	1.143	1.005	0.889	0.799	0.670	0.637	0.500	0.525	0.508	0.401	
438	0.414	0.395	0.423	0.417	0.423	0.447	0.471	0.479	0.465	0.508	
448	0.518	0.515	0.544	0.499	0.441	0.420	0.384	0.318	0.321	0.289	
458	0.265	0.227	0.193	0.156	0.150	0.142	0.130	0.122	0.118	0.114	
468	0.095	0.110	0.116	0.093	0.102	0.107	0.111	0.095	0.119	0.118	
478	0.109	0.098	0.088	0.089	0.088	0.079	0.081	0.090	0.068	0.050	_

Figure 7.9: Example of the text representation of a plot (emission scan of Anthracene).

7.4 As Contour Plot

Time resolved emission and transient absorption can be conveniently demonstrated as contour graphics as shown below. Like the 3D plot option, the Contour plot option is only available if the window contains more than one scan and all of them have equal step size (i.e. is a map).

The contour graphics with inactive cursor and active cursor are distinctively different. If the cursor is not active the whole window contains only the contour plot. On the other hand, if the cursor is active, two other data displays are shown beside the main contour graphic: a view from the x-axes and a view from the y-axes (see figure below). The axes view show the data of the active cursor position. Data can be highlighted by pressing the **Ctrl** button and the left mouse button at the same time. Highlighted data can be removed by holding the **Ctrl** button and pressing the left mouse button at the point of the intersection of the cursor cross hair.

The data in the axes windows can be normalised to the maximum by pressing the "N" button in the tool bar.

When the zoom function is chosen no cross hair is shown initially.

Additional contour lines can be added, contour lines can be removed, colours of contour lines can be changed, the resolution of the contour lines can be changed. Ref. to the contour plot options menu.



Figure 7.10: Example of a contour plot representation with active cursor (cross hair) and various marks

Contour plots options: The Contour plot options windows permit the selection of the colour properties for the background, axis and grid. In this type of plot, each of the experimental data is defined as an individual line with its own properties. In this way, you can add and remove specific lines from the plot just by clicking **Add** or **Remove** and modify, individually, the line settings and line width of each of them.

Current Contour Plot Opti	ons 🛛 🗙
Contour colour and line styles	Miscellaneous Options
Background Axis Grid Contour Line 1 Contour Line 2 Contour Line 3 Contour Line 4 Contour Line 5 Contour Line 6 Contour Line 7	Show Cursor Show Grid Normalise Cross Sections Scale Data From Zero Resolution 1: 1
Add <u>H</u> emove	
Contours = 68	
Line Width 1	Leset to Deraults Save as Defaults <u>Dk</u> <u>Cancel</u>

Figure 7.11: Contour plot options.

Miscellaneous options

Show cursor: This will enabled a Cursor in the window to read X and Y positions.

Show Grid: This will create a window grid.

Normalise Cross Sections: As default, if this option is not ticked, the Y scale for the side windows on the bottom and right of the main contour window will have a constant interval, where the maximum Y value is defined by the biggest trace. On the other hand, if the "Normalise Cross Section" option is ticked, the Y scale range will be re-scaled to the maximum Y-value for each curve. The same effect is obtained if the normalised button on the toolbar is clicked:

Resolution: Changes the resolution on the x-axes. This option is particular useful for decay data with large number of data channels. The editable value will define how many data points are averaged into one data point.

Scale Data from zero: As in 2D plot options.

Change Font: This will launch a "Font Dialog" box.

Reset to defaults: This will load the last default settings saved.

Save as defaults: This will save the current settings as default.

8 Data Analysis

The data menu is designed for processing the data acquired with the LP920. This menu has some common options for all the data type and others that are specific for spectral and lifetime scans. The menu will be only enabled if a measurement window is opened or if data has been acquired.

The general procedure is:



The following sections describe how to use the common and specific data manipulation options. The subtitles will indicate whether the data option is applicable for spectral data only, for lifetime data only, or whether they are applicable for both.

8.1 Combine (Lifetime and Spectral Data)

This option adds, subtracts, multiplies or divides scans.

- 1. Ensure the measurement window is in focus and contains only the scans that you want to work with. The operation will apply to all scans within the window so it may be necessary to extract a single scan. See next section about how to work with measurement windows.
- 2. On the **Data** menu, select **Combine**, and then click the required function.
- 3. On the list in the dialog box, select the scan to be added, subtracted, etc., from the list of all currently loaded scans.
- 4. If the measurement is not a map, choose whether you want the resultant scan(s) to be added to the same window or placed in a new window.

As an example, consider the addition of two traces called Sample1 and Sample2, where the measurement focus window is Sample2.

To add both traces together, the user has to click in the **Data menu/Combine/Add**. A dialog box will launch as shown in Figure 8.1.

The example in the dialogue box shows that Sample1 will be added to Sample2. The result can be shown in a new window **Create new window** or **Add to existing window**.

Add:	Measure	L900 Type	Range		
	Sample1	Time Scan	0.00 to 3999.60 ns		- 6
	🔺 Sample2	Time Scan	0.00 to 3999.60 ns		

Figure 8.1: "Add Data" dialogue box

A further function for combining data in the **Combine** submenu not mentioned above is **Append**. This feature allows the user to combine 2 plots of overlapping wavelength ranges. The function can only be applied to append one file to another, therefore it is only present in the file menu when the active window contains only one spectral scan. This function is especially useful when a spectral scan has been obtained in multiple parts (i.e. due to the use of second order filters). To append data, make sure the lower wavelength range scan is the active window and click **Data/Combine/Append**, next highlight the scan you wish the plot to be appended to. If there is an overlap of spectral ranges, the user can choose the wavelength at which the data is appended, by writing the value in the **Append at:** box (by default the software will insert the midpoint of the overlap range in this box). There may be an scaling offset (i.e. due to variation in transparency of different second order filters), this can be allowed for by clicking the '**Match amplitudes to**' tick box

Append Data					?×
Append	'Sample 3'				
	Contents of 'F90	10'			Ê
With:	Measurement	Туре	Range		
	Sample 3	Emission Scan Multiple Scans	365.00 to 406.00 365.00 to 500.00		
	🔄 Sample 4	Emission Scan	398.00 to 500.00		
	No overlap				
	nm	Г	Match amplitudes		
Results					
C Add to ex	kisting window		Create new	window	
			OK	Cancel	Apply

Figure 8.2: Append Data Dialog Box

Figure 8.3 shows an example of how to append two plots with overlapping spectral ranges and a mismatch in amplitude.



Figure 8.3: Procedure for appending plots.

8.2 Scale (Lifetime and Spectral Data)

This option allows re-scaling the data by a factor. The user has to click the measurement window that is going to be re-scaled (to make it active) and then go in the **Data/Combine/Scale** menu. A dialog box (Figure 8.4) will be launched, where the user can type the factor for re-scaling. The result can be displayed in a new window or an existing one.

S	Scale Data ? 🗙
	Carle (Carrelet)
	Scale Sample I
	by a factor of 0.5
	Results C Add to existing window © Create new window
	OK Cancel Apply

Figure 8.4: Scale Data Dialog Box

8.3 Normalise (Lifetime and Spectral Data)

ormalis	e Data to) Peak V	'alue			? >
Normalise p	eak of 'MultiS	icans'				
to	0.153968751	_				
Get Value f	rom Peak of:					
	Contents of '	Multi Scans'				- E
	Measure	Туре	Range			
	Sample2	Time Scan	0.00 to 3999.60 ns			
	🔺 Sample1	Time Scan	0.00 to 3999.60 ns			
	[⁻					
	J					
Baculto						
C Add to er	kisting window		Create	new win	dow	
			0	к	Cancel	Apply

Figure 8.5: Normalise Data Dialog Box

In order to normalise a scan(s), go to the option **Normalise** in the **Data** menu. This opens the **Normalise Data to Peak Value** dialog box (Figure 8.5). This option allows a value to be set for the highest peak in the spectral

scan, the rest of the scan is then scaled to fit. To do this type the value for the highest peak in the dialog box, set whether to add resultant plot to the current window or create a new window and click **OK** or **Apply**. In order to normalise the scan(s) to match the peak value of another scan, highlight it on the list in the dialog box, and its peak value will appear in the box. If you select a measurement container the maximum peak value from all its scans will appear, alternately, double-clicking on the measurement container folder will allow selection of a particular scan within the window.

Figure 8.6 shows an example of how to make the highest peak of 2 different plots the same height.





8.4 Subtract Baseline (Lifetime and Spectral Data)

This option allows a baseline to be subtracted from a plot. This is especially useful when using detectors with a high dark counts (e.g. near-infrared photomultiplier tubes). Click on **Data/Combine/subtract baseline** and in the dialog box (Figure 8.7), type the Y value to be subtracted from each point on scan. The resultant plot can be shown in a new window or added to an existing window by clicking the relevant radio button and clicking **Apply** or **OK**.

Subtract Base Line	? ×
From 'Sample 1'	
Subtract Line: Y = 10000	
Results C Add to existing window © Create new windo	w
OK Cancel A	۱pply

Figure 8.7: Subtract Baseline Dialog Box

8.5 Crop Range (Lifetime and Spectral Data)

Crop Data Range	?×
Crop the Range of 'Sample 2' To: 331.00 to 579.00	
Results O Add to existing window O Create new window	, ,
OK Cancel Ap	ply

Figure 8.8: Crop Data Range Dialog Box

This option allows the user to crop the X-range data from the active measurement window. Click on the **Data menu/Crop Range** and set up the maximum and minimum limits in the dialog box (Figure 8.8). When the crop range dialog box is opened, the current viewable wavelength range is entered into the crop range boxes.

Therefore, a easy way to crop the range of a scan is to zoom into the area of the scan that you want, press **Crop Range** then click **OK**. The x-axis will then be cropped to that in the magnified area.

8.6 Smooth (Spectral Data)

It allows smoothing the data by applying a binomial approximation. To smooth a plot, open the **Smooth** dialog box (Figure 8.9), which is accessed through the submenu **Data/Smooth**. Select the number of adjacent points that you want to take into account when smoothing (3, 5, 7 or 9) and click **OK**. To smooth with a larger number of points repeat process, e.g. smoothing again with 3 points will effectively add another two to the previous number of points.



Figure 8.9: Smooth Data Dialog Box

The smoothing routine used by the software is effectively a weighted average. For smoothing by 3 points, the software multiplies each point by 2, adds adjacent points and divides by 4 (the outermost points are multiplied by 3, added to there adjacent point and divided by 4). Smoothing by 5 points repeats this twice, smoothing by 7 points repeats it 3 times and smoothing by 9 points repeats it 4 times; This has the same effect as a weighted average of 5, 7 and 9 points.

8.7 Differentiate (Spectral Data)

This function calculates the derivative of n order of a active data scan. To perform a differentiation on a plot, open the **Differentiate** dialog box (Figure 8.10), which is accessed through the submenu **Data/Differentiate**. The function allows the user to differentiate from the 1st to the 9th order. If you want the scan(s) smoothed before differentiation, tick the **Smooth First** check box and select the number of points (see Smoothing, Section 8.6).

Calculating the derivative of a spectral scan computes the rate of change of the counts with respect to the wavelength. Figure 8.11 shows an example of a differential of an emission scan. Calculating n derivative orders is analogous to carrying out the differentiation n times, for example, differentiating a spectral scan to the second order finds the rate of change of the rate of change of the counts with respect to the wavelength.

Differentiate Data	? ×
Differentiate 'Sample 1' Smooth First Number of Points: 3 Order: 1st	
Add to existing window Create new window	

Figure 8.10: Differentiate Data Dialog Box



Figure 8.11: Spectral Scan and its Derivative

8.8 Integrate

This function calculates the integral of an active data scan. To perform an integration on a plot, open the **Integrate** dialog box (Figure 8.12), which is accessed through the submenu **Data/Integrate**. An integral function is used to find the area under a curve, the software creates a plot of the integral function by making each point equal to the integral of the plot to the left of it. Therefore, in order to find the area under the whole plot, the cursor (see Section 9.3.6) has to be used to find the integral value of the last point. Figure 8.13 shows the integral function of a spectral scan.

Integrate Data	? ×
Integrate 'Sample 1'	
Results O Add to existing window O Create new window	
OK Cancel Apply	,

Figure 8.12: Integrate Data Dialog Box



Figure 8.13 (a) Spectral scan, (b) Integral function

8.9 Reverse

This option performs the task of reversing the order of the data. It has the effect of producing a mirror image of the source data with the axis of symmetry being the midpoint of the data. The x-axis is not rescaled. This operation can be used to assist in comparisons of emission and absorption spectra. The dialog box used to carry out this operation is shown in Figure 8.14.

Reverse Data	? ×
➡ Reverse 'EmScan1' Data on X Axis	
C Add to existing window C Create new window	
OK Cancel App	ily

Figure 8.14: Reverse data edit box.

8.10 Peak Search (Spectral Data)

This option is designed to search for maxima and minima over a scan. To view the values of the peaks on a plot, click **View/Peaks**, or alternately click the icon on the toolbar:

In order to configure the 'show peaks' option to find peaks of the correct height, the **Peak Parameters** dialog box has to be accessed (Figure 8.15) through the submenu **Data/Peak Search Setup**.

The peak searching algorithm smoothes the data before searching for peaks and troughs. This technique helps to reduce the chance of noise values being wrongly detected as a peak feature. In the dialog box (Figure 8.15), the **Number of Points** parameter defines the level of smoothing. Peak features in the smoothed data set are discriminated by the **Shoulder Y Level**. This level details the minimum intensity for a smoothed feature to be recognised as a peak or trough. Features below this value will be ignored.

Peak Parameters	? ×
Standard Setting	Display
Shoulder Y Level: .10	🗖 Peak 🕜 Y Axis
Number of Points:	I Trough ⊙ XAxis
Additional Setting	
Sensitivity: 0	
Set Default OK	Cancel Apply

Figure 8.15: Peak Parameters Dialog Box

Either (or both) peaks and troughs can be marked in the measurement window. The mark used is a vertical line with either the x-axis position or the number of counts (y-axis) of the feature. Figure 8.16 shows a scan with peaks shown on it. The parameters were setup with the dialog box shown in Figure 8.15.

The **Sensitivity** value defines the specific noise amplitude to be used for peak rejection. Any feature with relative amplitude less than this value will be considered as noise. Note that the relative amplitude of a feature is its height above a baseline between the base points of the feature.



8.11 Correction (Spectral Data)

Once a spectral measurement has been acquired, if spectral correction was not carried out during acquisition, it can be corrected later on using the **Correction** option in the **Data** menu. Correction file measurements allow the spectral characteristics of the spectrometer to be measured so that scans can be adjusted to compensate for the transmission properties of the spectrometer at each step wavelength.

When there is a spectral file to be corrected, clicking the **Correction** option opens the **Manual Correction** dialog box (Figure 8.17).

1anual Correctio	n of Multi Scans : Multiple Scans
Correct sample 9 sample 8	Spectral Excitation Scan Correction
sample 7 sample 6 sample 5	Spectral Emission Scan Correction Emission Arm 1, Detector R928-P
· ·	Spectral Synchronous Scan Correction
Results	
C Add to existing	g window 💿 Create new window
	OK Cancel Apply

Figure 8.17: Manual Correction Dialog Box

9 Windows-related Menus

9.1 File Menu

The file menu makes available typical Windows facilities to let you open and save files, as well as import, export and print them. Also, some other features that are specific to Edinburgh Instruments software, such as Properties, are included in this menu.

File	Edit	View	Data	Options	Window I
M	lultiple	Measu	rements	5	
N	ew Sp	ectral M	leasure	ment	•
0	pen			Ctr	l+0
R	ecent	Files			•
S	ave			Ctr	l+s
S	ave As				
C	lose				
C	lose Al	I			
P	roperti	es			
Ir	nport f	rom Os	cillosco	pe Ctr	I+Alt+O
Ir	nport (SEM			
Ir	nport /	ASCII			
E	xport #	ASCII .			
P	rint Se	tup			
P	rint			Ctr	I+P
E	×it			Alt	+X

Figure 9.1: File menu layout.

New Spectral Measurement: The details of the measurements that can be performed with the F900 software are discussed fully in Chapter 5 (spectral) and Chapter 6 (lifetime).

Open: Clicking this menu item launches the **Open** dialog box that allows the user to select the data file name and then open it. In the F900 software, the data files have extension **.fl** for kinetic traces and **.fs** for spectral files. The dialog box is also accessible by clicking **File Open** icon in the toolbar.

Recent Files: This menu item allows the user to select one of the most recently opened files.

Save: This menu item allows the user to store the contents of an active data window under the current filename. A window becomes the active window simply by clicking on it. The data can be saved in one of two formats: EI measurement format or Windows Metafile format. The latter one can be easily copied and pasted into another applications such as word processors. It is also possible to have access to this menu by clicking the **Save** icon in the toolbar.

Save As: This launches a **Save As** dialog box to save the contents of the active window as a file whose name is created by the user. As with the save option, the data can be saved in one of two formats: EI measurement format or Windows Metafile format.

Close: This closes the active data window. The user will be prompted to save any unsaved data.

Close All: This closes all the windows displayed, no matter whether they are active, inactive, or minimised. Before any of them close, there will be a prompt to save them.

Properties: The scan properties are the set of measurement conditions of a measured scan. The F900 software can display the properties information, present them to be edited, and also to export them to printer, file and clipboard. The **Properties menu** includes a full description of all the conditions used to collect the data (time window, time calibration, etc). Most of them can be edited by clicking in the **Edit values** box. To display the scan properties on the file menu, click properties or double click the **Properties** icon.

Properties: The scan properties are a set of measurement conditions of a measured scan. F900 software can display the properties information (Figure 9.2), present them to be edited, and also to export them to printer, file and clipboard. The **Properties menu** includes a full description of all the conditions used to collect the data (time window, time calibration, etc) and most of them can be edited by clicking in the **Edit values** box. To display the scan properties perform one of the following options:

- On the File menu (or the right mouse click menu), click Properties.
- Double click the **Properties** icon.

If the measurement container has more than one scan, the screen shows information relevant to the container, and allows you to select one scan from **Measurement List** (see Figure 9.2). The properties of the scan are then displayed by pressing the **Properties** button, or by double clicking on the scan name in **Measurement List**. If only one scan is in the container, the scan properties' window is shown directly.

Timescan Container Properties ? 🗙
Description : Multi Time Scan
Type : Exponential Fit Time Scan
Comment :
Range (ns): 0.000 to 499.512
Measurement List:
Dil 1000 points
Oil 1000 pointsF1
Properties
OK Cancel

Figure 9.2: Example of Container Properties box (fit container)

The name of a scan, which has been selected from the measurement list, can be altered by typing. The alterations are shown 'live' on the graph. Pressing the **Cancel** button restores the original name(s). Scan properties can be edited at any time. The comment for the container file acts as the graph window's main title when the container is displayed.

Description : 🛛	Anthrascer	ne Emission	n			
Type : E	mission S	ican				
Comment :						
Correction Status : 🕅	No correct	tion				_
Em Wavelength :	365.00	to 500.0	0	step :	1.00	nm
Ex Wavelength :	357.00	nm				
Dwell time (s) :).20	per repeat		Repeats :	5	1
Lamp Source : 🕨	<e900< td=""><td></td><td></td><td>Temp(°C):</td><td>0.0</td><td>[</td></e900<>			Temp(°C):	0.0	[
-Scan Arm(Em) Par	ameters		Fixed	Arm(Ex) Par	ameters	
Polariser(*) : 🚺	No Polaris	er	F	olariser(*) :	No Polari	ser
Slit (mm) : 0).00	- 1		Slit [mm] :	0.00	_
Detector : F	1928-P					

Figure 9.3: Example of a scan properties page (spectral scan).

🛛 Impor	t ASCII file	s		? ×	
Convert files					
List of files	to import	"Sample data	".		
Type of co	Type of container to create		scans	•	
Data Delimit	er Character				
🔲 Tab	🔽 Comma	🔲 Space	🔲 Semicolon 🛛	Other: x	
🔽 Skip rep	eated delimiters				
Data starts	at row: 10 🍝	T Auto	o-detect for each file	•	
Line No.		1.0105.1	LOIS		
Row IU	0.00	-1.919E-1			
Row 12	0.200	-1.007E-1			
Bow 13	0.400	-1 900E-1			
Row 14	0.800	-1.880E-1		-	
Data Colum Import Dat	Data Columns Import Data from column: 1 🚔 🔽 Read X-values from column: 1 🛫				
		<u>0</u> K	<u>C</u> ancel	Apply	

Figure 9.4: Import ASCII dialog box

The contents of the window can be transferred to the clipboard by pressing the **To clipboard** button and printed by pressing the **Print** button. If the contents of the scan have been edited, they are checked before any action is taken. When using these functions the date of printing is attached. The format of the date string is the "Short Date style" specified on the **Date** tab of the Regional Settings icon in the Windows Control Panel.



Figure 9.5: ASCII format export dialog box

Import GEM: This option allows you to import data obtained with previous versions of the F900 software running under GEM operating system format (GEM is the data format that was used in spectrometers prior to 1995) into the F900 Windows version. The same effect is obtained by clicking the **Import Gem** icon.

Labels	Anthrascene Emission
Туре	Emission Scan
Comme	nt
start(nm) 365.00
stop(nm) 500.00
step(nm) 1.00
fixed/off	set(nm) 357.00
Xaxis	Wavelength
Yaxis	Counts
365.00	1.800E+3
366.00	2.806E+3
367.00	4.464E+3
368.00	7.520E+3
369.00	1.371E+4
370.00	2.214E+4
371.00	3.030E+4
372.00	4.269E+4
373.00	7.198E+4
374.00	1.335E+5

Figure 9.6: Section of data exported in ASCII format by F900 software

Import ASCII: This option allows the user to load data in ASCII format and perform a conversion into EI data format. Every time that a data file is open or imported from ASCII or GEM into L900, the software identifies if the data are kinetic traces, spectral scans or time-resolved data and creates a specific container according to this. A series of check boxes on the dialog menu (Figure 9.4) allows the user to choose the type of delimiter and which columns contain the data for each axis. This option can also be accessed by pressing the **Import ASCII** toolbar button.

Export ASCII: This option allows the data to be exported in ASCII format and subsequent importing of the data into other applications (such as spreadsheets) that use the ASCII format. When the **Export ASCII** option is selected in the File menu or when the **Export ASCII** icon is clicked, the system launches a **Data export format** box (Figure 9.5). A pair of radio buttons permit the user to choose whether the data is to be saved in rows or columns. In addition, a set of radio buttons allow the user to choose a character that will serve as a separator between the numeric values in the raw data. The importing application can then be configured to recognise this separator (or delimiter) and, in the case of a spreadsheet, display the data in a suitable configuration of rows and columns. The properties can also be included in the exported file of the measurement by clicking in the **Include properties** box. These properties will be located at the beginning of the ASCII file. An example of the exported data is shown in Figure 9.6.

Print setup: This launches a standard Print Setup window.

Print: This launches a Print Dialog Box as shown in the picture below. The user can select the colours to be used by the printer to plot the data (Printer Colour Options) or select to print it as it is shown in the screen (Use Screen Colours).

The print option is also accessible by clicking the **Print** icon.

Print	×
Printer: Panasonic KX-P6300 GDI	
Printer Colour Options	Miscellaneous
Background	Print Properties
	Print Annotation
Colour Change	Print Logo
Line Style Line, Solid	Print Title
	Print Grid
Line Width 1	
Setup Prin	nt Cancel

Figure 9.7: Print dialog box

Exit: This closes the F900 software. The user will be prompted to save all the data unsaved.

9.2 Edit Menu

The only option available in the Edit menu is a **Copy** (Ctrl-C) facility that allows copying of the active window, and (Ctrl-V) to paste it in another program.

9.3 View Menu

Figure 9.8: View menu options.

When the View Menu is selected, the drop-down box shown in Figure 9.8 is displayed. The various items in this menu will now be described.	µF Lamp Setup Signal Rate
	 As 2D Plot As Contour Plot As 3D Plot As Text
	Plot Options Zoom In Cursor Grid Peaks Normalise
	✓ Status Bar✓ Tool Bar

9.3.1 Lamp setup

This menu item is used to setup the lamps installed with the system. Depending on the configuration, either or both of the microsecond flashlamp and the nanosecond flashlamp may be available.

µF Lamp Setup	? ×
Frequency: 100.0	♣ Hz
OK	Cancel

Figure 9.9: Dialog box to set up the microsecond flashlamp.

The dialog box used to set up the microsecond flashlamp is shown in Figure 9.9. The only option that can be set here is the lamp frequency. This has a resolution of 0.1 Hz and can be set in the range 0.1 - 100.0 Hz.

The dialog box for the nanosecond flashlamp is shown in Figure 9.10. The settings that can be made here are closely interrelated and reference should be made to the "nF 900 Nanosecond Flashlamp" operating instructions manual for compatible values of EHT, Frequency, Gas and Gap. Pressing Switch Lamp On will start the lamp and the parameters in the Measured Parameters grouping will be activated. The Auto Start check box is used to indicated whether the nanosecond flashlamp should start as soon as a measurement sequence is commenced.

nF Lamp	? X
Set Up	1
EHT: 6.50 🛨 KV	₽
Frequency: 40.000 🚔 kHz	Switch Lamp Un
	Auto Start
Measured Parameters FHT (NA)	6 50 10
	0.00
0 Frequency (kHz)	0 250 1
	0.200
0 Pressure (bar)	0.00 2
Stored Parameters	
Gas: Hydrogen 💌	Gap: 1.00 🚔 mm
	Apply

Figure 9.10: Dialog box for setting up the nanosecond flashlamp.

9.3.2 Signal Rate

Selecting this menu item opens the Signal Rate dialog box that is described in Section 5.1.

9.3.3 As 2D Plot, As Contour, As 3D Plot, As Text

These options are used to set how the data is viewed and they are controlled by a radio button so that only one of them is valid at any particular time. The **As Contour** and **As 3D Plot** options are only available for map measurements. Further details of these settings are provided in Chapter 7. The currently selected plot style is indicated by a black dot next to it.

9.3.4 Plot options

Selection of this menu item open the **Plot Options** dialog box that is used to control how the data is presented. Further detail of plot options are described in Chapter 7.

9.3.5 Zoom In

Selecting this menu item places a cross-hair cursor on the plot. Dragging this cross-hair will produce a rectangle and when the mouse button is released the plot will zoom-in to the area delineated by this rectangle. Once a plot has been zoomed, a **Zoom Out** menu item will also appear. Selecting this menu item will restore the plot to its original scaling. If a plot is already zoomed the nan arrow is displayed next to the menu item.

9.3.6 Cursor

Selecting this option will place a vertical cursor on the plot. Moving this cursor with the mouse will cause the co-ordinates of the data point beneath the cursor to be displayed. The x-ordinate is displayed at the bottom left of the plot and the y-ordinate(s) is/are displayed on the right hand side of the plot. The menu item has a tick mark next to it if the cursor is visible.

9.3.7 Grid

Selecting this option will superimpose a grid on the plot. The presence of a grid is denoted by a tick mark next to the menu item.

9.3.8 Peaks

Selection of this item will display the amplitudes of the peaks that are present in the plot. The criteria for peak selection are made with the **Peak Search Set Up** menu item described in Section 8.10.

9.3.9 Normalise

This option is used with contour plots to normalise the data to the selected data value. The use of this option is indicated by a tick mark next to the menu item.

9.3.10 Status bar

Selecting this option will display the status bar that shows the current settings for the spectrometer hardware and also messages that have been reported by the system. This visibility of the status bar is indicated by a tick mark next to the menu item.

9.3.11 Tool bar

Selecting this option will display the toolbar containing the Edinburgh Instruments speed buttons. The visibility of the toolbar is denoted by a tick mark next to the menu item.

9.4 Options Menu

When the **Options Menu** is selected then the drop-down menu shown in Figure 9.11. These options will now be discussed in more detail.

Plot Defaults...
Correction Files...
Window Options...

Hardware Configuration... PCS900 Configuration... Drivers & Cards...

Figure 9.11: Options menu items.

9.4.1 Plot defaults

This menu item opens a dialog box containing the plot settings for all the measurement types available for the F900 software. Further details of plot settings are provided in Chapter 7.

9.4.2 Correction files

If it is intended to correct captured data for the effects of instrumentation then correction files have to be used. To inform the system of the files to be used, click **Correction Files**. The dialog box shown in Figure 9.12 will then be displayed.

Correction Files	? ×
Ref. Detector Excitation Emission1	
Reference Detector File	
H:\Test Data\Correction Files\Test-correction.FS	Modify
Range 200.00nm to 900.00nm, step 1.00nm	
ОК	Cancel

Figure 9.12: Correction files selection dialog box.

This dialog box permits correction files to be entered for the Reference Detector, the Excitation monochromator and one or more Emission monochromators. The tab for the Reference detector is shown in Figure 9.12, but the other two tabs are identical. To enter a file name press the **Modify** button. This will open a Windows **File Open** dialog box. Navigate to the correct directory and file and then press **Open**. This will place the file name in the **Reference Detector File** edit box. If no correction file is to be applied to a measurement then the **Reference Detector File** edit box should be cleared.

9.4.3 Windows options

The **Windows Options** dialog box is used to select the methods of handling the plot windows. It is accessed by pressing the **Windows Options** item in the **Options** menu.

Window Options 🛛 🔋 🗙	
 Join multiple files on open Append filenames to scans on combining Close source window on split Close source window on extract Close source windows on join visible Autocorrect old files on open Recently used file list: 5 = entries. Temperature Tolerance +/- 0.00 = °C/K 	
OK Cancel Apply	

Figure 9.13: Dialog box for windows options.
The Items that can be checked are:

Join multiple files on open: If this option is selected then when multiple files are selected with the **Open** menu (Section 9.1) they will, if possible, be loaded into a single plot. Otherwise they will be loaded into individual plots. Files can only be loaded in to one plot if the are the same measurement type.

Append filenames to scans on combining: If this option is selected then previously saved files will have their filename appended to their scan name when they are involved in combine operations.

Close source window on split: If this option has been selected then the source window for the waveforms will be closed when a split operation is performed, so that only the newly created windows will be displayed.

Close source window on extract: If this option has been selected then the source window for the waveforms will be closed when an extract operation is performed, so that only the newly created windows will be displayed.

Close source windows on join visible: If this option has been selected then the source window for the waveforms will be closed when a join visible operation is performed, so that only the newly created windows will be displayed.

Autocorrect old files on open:

The other items in this dialog box are:

Recently used file list: The number of files to be explicitly displayed in the File drop-down menu.

Temperature Tolerance: When separate files are combined, checks are performed to see whether the combined files constitute a map measurement. One of the checks is on whether the temperatures are the same. This check compares exact temperature values so a small difference will prevent a map being established. Entering a non-zero value in this edit box will make the comparison process examine whether there is an overlap in the temperatures when the tolerance value is taken into account. If so, then a map will be allowed.

9.4.4 Hardware configuration

This menu item opens the Hardware Configuration dialog box that is described in Chapter 3.

9.4.5 PCS900 configuration

PCS900 Configuration	×
Threshold DACs	
Ch 0: <u>19.7</u> 🌩 mV Ch 3: 19.7 🌩 mV	
Ch 1: -19.7 🔹 mV Ch 4: 250.0 🔹 mV	
Ch 2: -19.7 🔺 mV Trig: 75.4 🔺 mV	
Version Information	
DLL Version V2.0	Οκ
Hardware Version V1.0	
Hardware Type PCI Bus	Cancel

Figure 9.14: PCS900 configuration dialog box.

Selecting the **PCS900 Configuration** option will open the dialog box shown in Figure 9.14. This box permits changes to be made to the DAC (Digital to Analogue Converter) thresholds on each of the card channels. In addition, software and hardware version information is supplied for the user.

9.4.6 Drivers and cards

Selecting this option will open an 'information-only' dialog box, shown in Figure 9.15. The information contained in the box will show the presence of the DLLs used to interface to particular hardware component and also the presence of the hardware component.

Configuration Summary					×		
	Detected Elements						
	V	TCC-900 DLL	$\overline{ } $	TCC-900 Card			
	Г	Norland DLL	Г	Norland Card			
	$\overline{\mathbf{v}}$	PCS-900 DLL	$\overline{ \mathcal{V} }$	PCS-900 Card			
	$\overline{\mathbf{v}}$	CD900 DLL					
	$\overline{\mathbf{v}}$	Level1 DLL					

Figure 9.15: Summary of the presence of hardware options and of the software utilities used to interface to them.

9.5 Window Menu

The window menu is divided in three parts: the first one is a general feature to all Windows environments and includes options to display the windows vertically, horizontally or in cascade.

Tile Vertically: to fit all open windows into the client area with their heights maximised.

Tile Horizontally: to fit all open windows into the client area with their widths maximised.

Cascade: to overlap all open windows with their title remaining visible.

Arrange lcons: to arrange all minimised windows so that their icons don't overlap.

When tiling or cascading make sure all the windows to be displayed are open. Minimised windows cannot be displayed.

The second part of the Window menu is specific for the F900 software and includes options to display and work with data windows.

Note: Every time that the user wants to do some operation with a data window, it is first necessary to set up the focus of input in that window. To do this, perform one of the following two procedures:

- select the name of the window to be made active in the list.
- left click the mouse on any part of the window.

The operations that can be performed on the data windows are:

Extract: This operation extracts one or more scans from a window that contains more than one scan. To perform the operation follow these steps: in the **Window** menu, click **Extract** and select the scans to be extracted from the measurement list. Alternatively, right click in the active Window and select **Extract**. A further alternative for running this operation is to press the **Extract** toolbar button.

Split: The Split function splits a window with more then one data set into individual scans. To perform the operation go to the **Window** menu and click **Split**. Each scan in the source container will be displayed in its own window. Alternatively, right click in the active Window and select **Split**. A further alternative for running this operation is to press the **Split** toolbar button.

Join Visible: This option joins separate data windows into a single one, provided the data are of the same type, e.g. windows containing spectral data and other windows containing kinetic data can not be joined together. The function does not have any effect on minimised data windows. On the **Window** menu click **Join**. Compatible scans in all the open windows will be merged into a new window. An alternative for running this operation is to press the **Join Visible** toolbar button.

Insert scan: This option allows the addition of a scan into a data window. After selecting this function a list of all windows and the individual data within those windows is provided. From this, the scans to be inserted can be chosen. To perform this operation, select **Insert scans** from within the **Window** menu. This function is also available by right clicking in the window and selecting **Insert scans**.

Delete Scan: This option permits the deletion of a specific scan from a data window. Deleted data will not be retrievable. To perform the operation, click **Delete scans** in the **Window** menu. Alternatively, right click in the Window and select **Delete scans**.



Figure 9.16: Graphical demonstration of some important Window manipulating functions.

Additional comments about working with data windows.

During the operation of fluorescence spectrometers, many data sets will be created, and manipulation of these data sets can generate a large number of data windows. In order not to get lost among the many data windows it is advisable to minimise as many data windows as possible and to title the windows appropriately.

9.6 Help Menu

The help menu contains two items, which are

Help Topics: Pressing this menu item opens a "windows style" help dialog box containing information that is designed to assist the user in using the F900 software.

About F900: This dialog box displays version information regarding the currently installed F900 software.

10Warranty

- 1 a) The Company guarantees the equipment forming the subject of the contract/quotation against defective materials and workmanship for a period of one year from the date of delivery to the Purchaser.
 - b) In the case of sub-assemblies of equipment not manufactured by the Company, but incorporated in the equipment ordered, the Purchaser will be entitled only to the benefit and/or limitations of any guarantee given by the makers of such assemblies.
 - c) In no event shall the Company be liable for any consequential loss or damage arising from failure of the equipment under warranty.
 - d) At the end of the one year period referred to herein, all claims upon all liability of the Company shall be absolutely at an end.
- 2 a) The Company also warrants that the equipment conforms to specifications confirmed in writing at the time of order acknowledgement.
 - b) No warranty is made or implied as to the suitability of any equipment for the Purchaser's intended use beyond such performance specifications as form part of the contract.
- 3. The purchaser warrants:
 - a) That he will carefully examine and list all parts of the equipment supplied by the Company and notify the Company in writing of any shortage, defect or failure to comply with the contract, which is or ought to be apparent upon such examination and test, within 48 hours of the equipment being delivered to or collected by the Purchaser.
 - b) The equipment will be operated in accordance with the instructions and advice detailed in the appropriate operating instructions manual, or any other instructions which may be provided by the Company. The Company shall not be held responsible for any defect arising from the Purchaser's failure to comply with these recommendations and instructions or from damage arising from negligence or exposure to adverse environmental conditions.
- 4. The warranty is effective when:
 - a) Any defects in the equipment supplied are notified immediately by the Purchaser to the Company.
 - b) The equipment is returned to the Company at its Edinburgh premises, transportation and insurance prepaid, and undamaged by the failure to provide sufficient packaging.
 - c) The Purchaser has made payment in full for the contract in accordance with the Company's normal trading terms, i.e. 30 days from date of invoice.
- 5. The warranty covers:
 - a) Engineer's time costs during inspection and repair.
 - b) Any materials or components, which require to be replaced.
 - c) Return carriage costs to the Purchaser
- 6. However, if the Purchaser requests a service engineer to carry out the necessary inspection and repair of the equipment covered by the warranty on site, the Purchaser will be liable, at the Company's discretion, for:
 - a) Engineer's travelling time costs.
 - b) Engineer's travelling and accommodation expenses.

The timing of the inspection and repair of the equipment will be determined entirely at the discretion of the Company.

11CE Declaration of Conformity

CE

Manufactured by:

Edinburgh Instruments Ltd. 2 Bain Square Livingston EH54 7DQ Tel.: + 44 1506 425 300 Fax.: + 44 1506 425 320

Applicable Standards:

Generic Immunity Generic Emission Electrical Safety Standards EN 50082-1 : 1992 EN 50081-1 : 1992 EN 61010-1 : 1993

Edinburgh Instruments Ltd. certify that this equipment conforms to the protection requirements of the above Directives.